E Instructions for the interior body exchange of the CO₂ sensors InPro® 5000 (i)

This instruction sheet explains how to replace the interior body (pH electrode) of the InPro 5000 (i) sensor. The information on this sheet and the safety requirements in particular, must be observed by the operator and all other persons working with this equipment.

Order no. of replacement packages

	for InPro 5000	for InPro 5000 i
Interior body 120	52 206 057	30 019 049
Interior body 220	52 206 084	30 019 170
Interior body 320	52 206 085	30 019 175

Important notes:

- ☞ Great care should be exercised when handling the interior body which is made partly of glass, as any cracks will adversely affect sensor performance.
- ☞ Due to differences in sensor construction, the interior bodies of the InPro 5000 and InPro 5000 i cannot be exchanged.

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E The replacement package of the interior body contains:

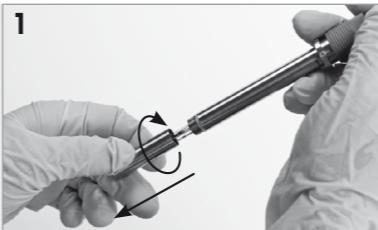
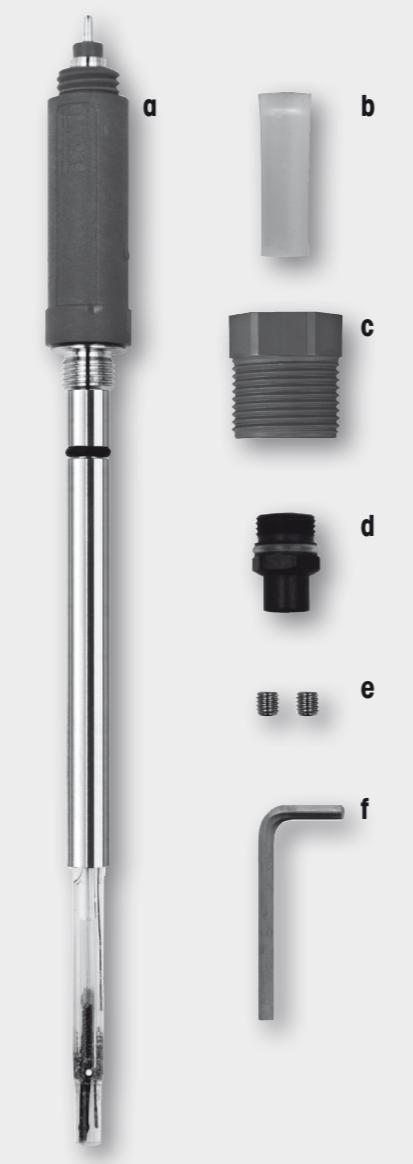
- 1 interior body (pH electrode) of specified length (a)
- 1 wetting cap (b)
- 1 Pg 13.5 threaded sleeve (c)
- 1 retainer nut (d)
- 2 hex screws (e)
- 1 allen wrench (f)

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E Instructions for the interior body exchange of the CO₂ sensors InPro® 5000 i

- 1 Unscrew the cap sleeve from the shaft and carefully pull it from the sensor.
- 2 Remove the membrane body. Usually the membrane body remains in the cap sleeve and should be ejected by pushing it with a flat finger tip. Otherwise pull the membrane body from the interior body. Rinse glass electrode with deionized water.

- 3 Unlock the interior body by unscrewing the hex screw counter clockwise with an allen wrench.
- 4 Remove the interior body by turning the upper part of interior body counter clockwise. Pull the interior body from the steel shaft. The used interior body, the Pg 13.5 threaded sleeve (red), the retainer nut and the hex screw cannot be used further. Please discard according to local waste regulations.

- 5 Mount the new Pg 13.5 threaded sleeve and the retainer nut to the steel shaft.
- 6 Insert the new interior body into the steel shaft. Turn clockwise and tighten until the interior body is fully engaged in the steel shaft.
- 7 Insert new hex screw and lock interior body with the allen wrench. Verify that the interior body is correctly engaged. For the next steps e.g. calibration, refer to the instruction manual of the InPro 5000 i CO₂ sensor.

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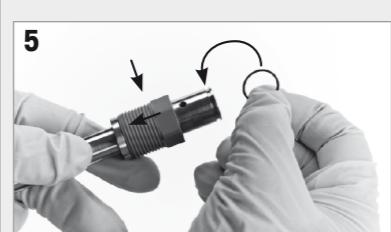
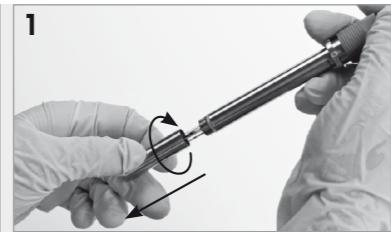
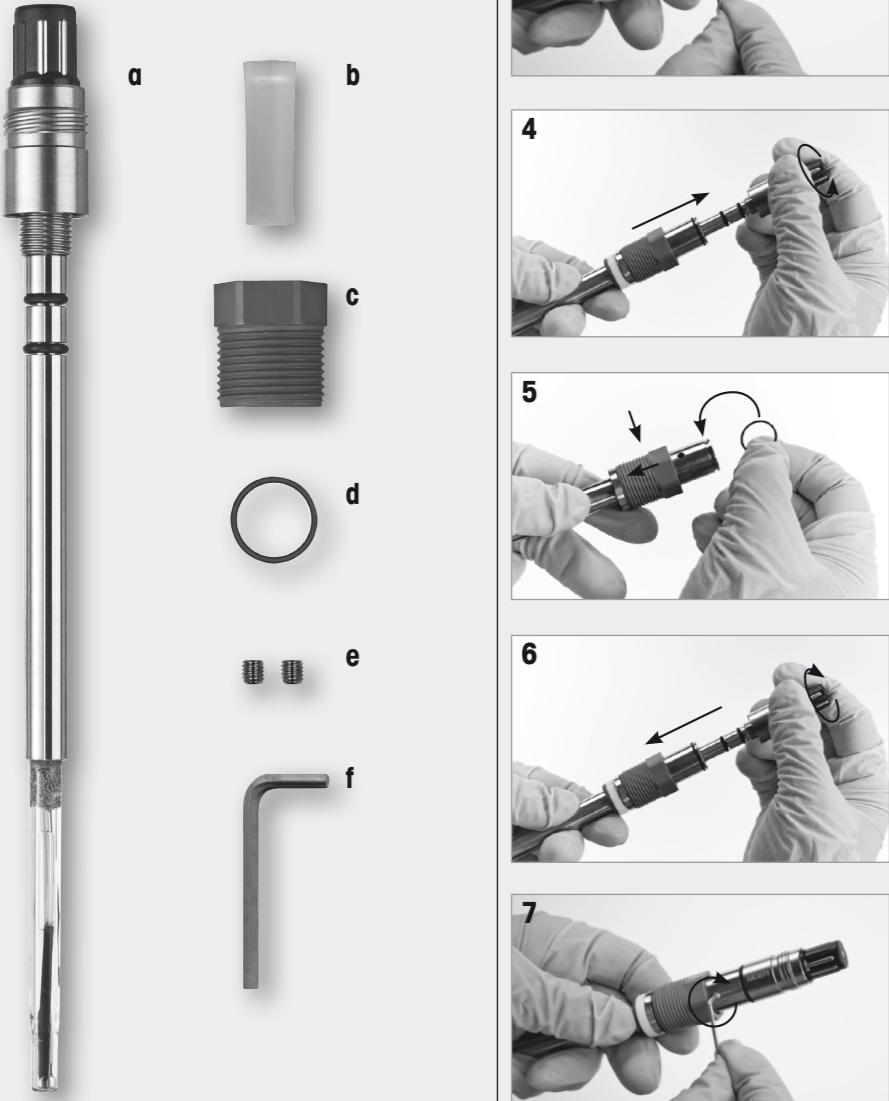
E The replacement package of the interior body contains:
 1 interior body (pH electrode) of specified length (**a**)
 1 wetting cap (**b**)
 1 Pg 13.5 threaded sleeve (**c**)
 1 O-ring, 13.0 x 1.0 mm, Viton® [0.51" x 0.04"] (**d**)
 2 hex screws (**e**)
 1 allen wrench (**f**)

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E Instructions for the interior body exchange of the CO₂ sensors InPro® 5000

- 1 Unscrew the cap sleeve from the shaft and carefully pull it from the sensor.
- 2 Remove the membrane body. Usually the membrane body remains in the cap sleeve and should be ejected by pushing it with a flat finger tip. Otherwise pull the membrane body from the interior body. Rinse glass electrode with deionized water.
- 3 Unlock the interior body by unscrewing the hex screw counter clockwise with an allen wrench.
- 4 Remove the interior body by turning the upper part of interior body counter clockwise. Pull the interior body from the steel shaft. The used interior body, the Pg 13.5 threaded sleeve (red), the O-ring and the hex screw cannot be used further. Please discard according to local waste regulations.
- 5 Mount the new Pg 13.5 threaded sleeve and the O-ring to the steel shaft.
- 6 Insert the new interior body into the steel shaft. Turn clockwise and tighten until the interior body is fully engaged in the steel shaft.
- 7 Insert new hex screw and lock interior body with the allen wrench. For the next steps e.g. calibration, refer to the instruction manual of the InPro 5000 CO₂ sensor.

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Analog Sensor InPro 5000

52 002 466

Mettler-Toledo AG, Process Analytics, Im Hackacker 15, CH-8902 Urdorf, Switzerland, Tel. +41 44 729 62 11, Fax +41 44 729 66 36
 Subject to technical changes. 10/11 © Mettler-Toledo AG. Printed in Switzerland

BR Mettler-Toledo Ind. e Com. Ltda., Alameda Araguado, 451 – Alphaville, BR – 06456-000 Barueri/SP, Phone +55 11 4166 74 44, Fax +55 11 4166 74 01

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Instructions for interior body exchange of the CO₂ sensors InPro® 5000 (i)

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D Mettler-Toledo GmbH, Prozeßanalytik, Ockerweg 3, D – 35396 Giessen, Phone +49 641 507 333, Fax +49 641 507 397