
English	Reference Manual QLX3 Liquid Module for XPR Analytical Balances
Deutsch	Referenzhandbuch QLX3-Flüssigdosiermodul für XPR-Analysenwaagen
Español	Manual de referencia Módulo de líquido QLX3 para balanzas analíticas XPR
Français	Manuel de référence Module pour liquides QLX3 pour balances d'analyse XPR
简体中文	参考手册 QLX3液体模块 适用于XPR分析天平
日本語	リファレンスマニュアル QLX3液体モジュール XPR分析天びん用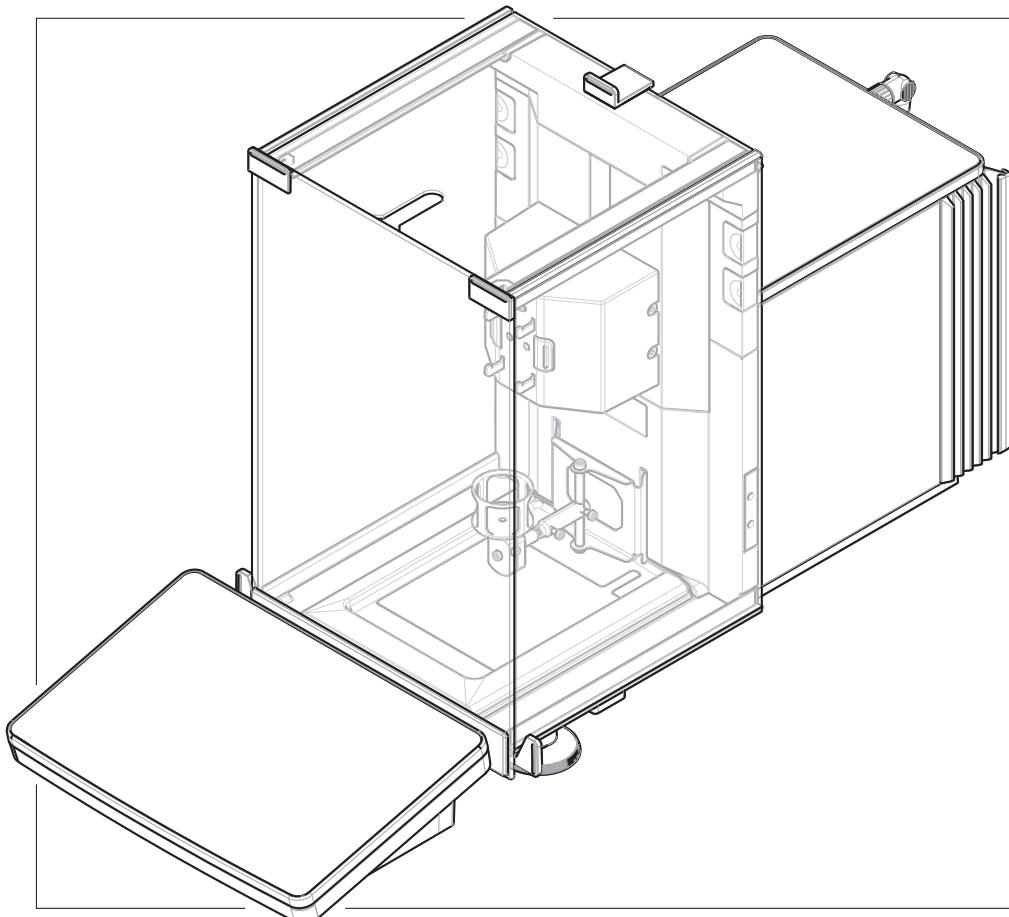

METTLER TOLEDO

Reference Manual **QLX3 Liquid Module**

English

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Table of Contents

1	Introduction	3
1.1	Further documents and information	3
1.2	Explanation of conventions and symbols used	3
1.3	Acronyms and Abbreviations	4
1.4	Compliance information.....	4
2	Safety Information	6
2.1	Definitions of signal words and warning symbols	6
2.2	Product-specific safety notes	6
3	Design and Function	10
3.1	Function description	10
3.2	Overview.....	10
3.3	Components description	10
3.4	Dosing head RFID tag	12
4	Installation and Putting into Operation	13
4.1	Scope of delivery	13
4.2	Selecting the location	13
4.3	Setting up the liquid dosing system.....	14
4.3.1	Interface of the pump	14
4.3.2	Wiring the pump	14
4.3.3	Connecting the tubes	14
5	Operation	19
5.1	Liquid dosing	19
5.1.1	Installing and removing the liquid dosing head.....	19
5.1.2	Adjusting the dosing height	19
5.1.3	Attaching the dosing head to the bottle cap	20
5.1.4	Handling the bottle	20
5.1.5	Using the QL3 pump	21
5.2	Using other weighing methods	22
6	Maintenance	23
6.1	Cleaning	23
6.1.1	Cleaning the housing.....	23
6.1.2	Cleaning the weighing chamber	23
6.1.3	Cleaning the liquid dosing elements	23
6.2	Replacing the sealing ring and fastening nut on the liquid tube	24
7	Troubleshooting	26
7.1	Error symptoms	26
8	Technical Data	28
8.1	General data	28
8.2	Model-specific data.....	28
8.3	Material specifications	29
8.3.1	Glossary of material acronyms	29
8.3.2	Liquid dosing head QL001	30
8.3.3	Liquid dosing head QL003	31
8.3.4	QLL standard/advanced kit for bottles	31
8.4	Dimensions.....	32

9	Disposal	33
10	Accessories and Spare Parts	34
10.1	Accessories.....	34
10.2	Spare parts	38
10.2.1	QLX3 liquid module.....	38
10.2.2	QL3 pump.....	38
10.2.3	QLL spare parts kit for bottles	39
Index		41

1 Introduction

Thank you for choosing a METTLER TOLEDO instrument. The instrument combines high performance with ease of use.

1.1 Further documents and information

► www.mt.com/XPR-automatic

This document is available in other languages online.

► www.mt.com/QLX3-RM

Instructions for cleaning a balance: "8 Steps to a Clean Balance"

► www.mt.com/lab-cleaning-guide

Search for software downloads

► www.mt.com/labweighing-software-download

Search for documents

► www.mt.com/library

For further questions, please contact your authorized METTLER TOLEDO dealer or service representative.

► www.mt.com/contact

1.2 Explanation of conventions and symbols used

Conventions and symbols

Key and/or button designations and display texts are shown in graphic or bold text, e.g., Edit.

Note

For useful information about the product.

Refers to an external document.

Elements of instructions

In this manual, step-by-step instructions are presented as follows. The action steps are numbered and can contain prerequisites, intermediate results and results, as shown in the example. Sequences with less than two steps are not numbered.

■ Prerequisites that must be fulfilled before the individual steps can be executed.

1 Step 1

 ⇒ Intermediate result

2 Step 2

 ⇒ Result

1.3 Acronyms and Abbreviations

Original term	Explanation
EMC	Electromagnetic Compatibility
FCC	Federal Communications Commission
LPS	Limited Power Source
POM	Polyoxymethylene
RFID	Radio-frequency identification
RM	Reference Manual
sd	Standard deviation
SELV	Safety Extra Low Voltage
SOP	Standard Operating Procedure
UM	User Manual
USB	Universal Serial Bus

1.4 Compliance information

National approval documents, e.g., the FCC Supplier Declaration of Conformity, are available online and/or included in the packaging.

► <http://www.mt.com/ComplianceSearch>

Contact METTLER TOLEDO for questions about the country-specific compliance of your instrument.

► www.mt.com/contact

United States of America

This equipment has been tested and found to comply with the limits for a **Class A** digital device, pursuant to part 15 of the FCC Rules. These limits are designed to provide reasonable protection against harmful interference when the equipment is operated in a commercial environment. This equipment generates, uses, and can radiate radio frequency energy and, if not installed and used in accordance with the instruction manual, may cause harmful interference to radio communications. Operation of this equipment in a residential area is likely to cause harmful interference in which case the user will be required to correct the interference at his own expense.

This device complies with Part 15 of the FCC Rules. Operation is subject to the following two conditions:

1. This device may not cause harmful interference.
2. This device must accept any interference received, including interference that may cause undesired operation.


Changes or modifications not expressly approved by the party responsible for compliance could void the user's authority to operate the equipment.

Canada

This device contains licence-exempt transmitter(s)/receiver(s) that comply with Innovation, Science and Economic Development Canada's licence-exempt RSS(s). Operation is subject to the following two conditions:

1. This device may not cause interference.
2. This device must accept any interference, including interference that may cause undesired operation of the device.

Singapore

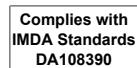South Africa

Thailand

2 Safety Information

Two documents named "User Manual" and "Reference Manual" are available for this instrument.

- The User Manual is printed and delivered with the instrument.
- The electronic Reference Manual contains a full description of the instrument and its use.
- Keep both documents for future reference.
- Include both documents if you transfer the instrument to other parties.

Only use the instrument according to the User Manual and the Reference Manual. If you do not use the instrument according to these documents or if the instrument is modified, the safety of the instrument may be impaired and Mettler-Toledo GmbH assumes no liability.

2.1 Definitions of signal words and warning symbols

Safety notes contain important information on safety issues. Ignoring the safety notes may lead to personal injury, damage to the instrument, malfunctions and false results. Safety notes are marked with the following signal words and warning symbols:

Signal words

DANGER A hazardous situation with high risk, resulting in death or severe injury if not avoided.

WARNING A hazardous situation with medium risk, possibly resulting in death or severe injury if not avoided.

CAUTION A hazardous situation with low risk, resulting in minor or moderate injury if not avoided.

NOTICE A hazardous situation with low risk, resulting in damage to the instrument, other material damage, malfunctions and erroneous results, or loss of data.

Warning symbols

General hazard

Notice

2.2 Product-specific safety notes

Intended use

This dosing system is designed to be used in analytical laboratories by trained staff. The dosing system is intended for weighing and dosing powder or liquid samples.

Any other type of use and operation beyond the limits of use stated by Mettler-Toledo GmbH without consent from Mettler-Toledo GmbH is considered as not intended.

Responsibilities of the instrument owner

The instrument owner is the person holding the legal title to the instrument and who uses the instrument or authorizes any person to use it, or the person who is deemed by law to be the operator of the instrument. The instrument owner is responsible for the safety of all users of the instrument and third parties.

Mettler-Toledo GmbH assumes that the instrument owner trains users to safely use the instrument in their workplace and deal with potential hazards. Mettler-Toledo GmbH assumes that the instrument owner provides the necessary protective gear.

Protective equipment

Chemical-resistant gloves

Goggles

Lab coat

Safety notes**⚠ WARNING****Death or serious injury due to electric shock**

Contact with parts that carry a live current can lead to death or injury.

- 1 Only use the METTLER TOLEDO power cable and AC/DC adapter designed for your instrument.
- 2 Connect the power cable to a grounded power outlet.
- 3 Keep all electrical cables and connections away from liquids and moisture.
- 4 Check the cables and the power plug for damage and replace them if damaged.

**⚠ WARNING****Injury and/or damage due to hazardous substances**

Chemical, biological or radioactive hazards can be associated with the substances processed by the instrument. During dosing procedures, small amounts of the dosed substance may become airborne and penetrate the instrument or contaminate its surroundings.

The substance characteristics and related hazards is the full responsibility of the instrument owner.

- 1 Be aware of possible hazards associated with the substance and take adequate safety measures, e.g., those stated on the safety data sheet provided by the manufacturer.
- 2 Make sure that every instrument part in contact with the substance will not get altered or damaged by the substance.

**⚠ WARNING****Injury or damage due to powder handling**

Powders may get compacted in the dosing head and block it. Too much force applied by the dosing head mechanism may lead to a break and potentially hazardous substances may get airborne.

- 1 Handle the dosing heads with care.
- 2 If the dosing head seems blocked, avoid further dosing. Remove the head from the instrument and turn it upside down to loosen the powder.
- 3 Cease work immediately in the event of leak or break.

**⚠ WARNING****Injury and/or damage due to reacting, flammable, or explosive substances**

During the dosing procedure, substances could be combined and cause an exothermic reaction or explosion. This includes powders, liquids, and gases.

The sample characteristics and related hazards is the full responsibility of the instrument owner.

- 1 Be aware of possible hazards associated with reacting, flammable, or explosive substances.
- 2 Ensure a working temperature low enough to prevent the formation of flames or an explosion.

⚠ WARNING

Injury or death due to toxic, explosive, or flammable substances

If you use toxic, explosive, or flammable liquids with the pump, the exhaust air will be contaminated.

- Connect a tube to the exhaust air outlet to collect the contaminated air.

⚠ WARNING

Injury and/or damage due to reacting substances

When pressure is released from the bottle, the air/gas in the bottle moves back towards the pump. The air/gas coming from the coupled outlets mixes in the pump. Molecules of the substances in the various bottles can get in contact through this contaminated air/gas.

- 1 Do not connect bottles with incompatible liquids to the same pump simultaneously.
- 2 Before connecting a second, incompatible liquid to the pump, disconnect the first bottle and purge the pump cavity with clean air/gas.

⚠ WARNING

Injury and damage to the pump or the bottle due to high pressure

High pressure from external gas can damage the pump or the bottle.

- 1 Use a regulator on the external gas line.
- 2 Ensure that the pressure of the external gas does not exceed 0.5 bar (7.2 psi).

⚠ CAUTION

Injury due to splashing liquids

If the pressure in the bottle is not released, liquid might splash when removing the micro dosing valve, opening the bottle, or removing the liquid tube.

- Always release pressure before removing the micro dosing valve, opening the bottle, or removing the liquid tube.

⚠ CAUTION

Injury due to leaking liquids

Wrongly cut tubing can result in leaking connections.

- Cut the tubes with a tube cutter or a sharp knife.

⚠ CAUTION

Injury due to moving parts

- Do not reach into the working area while parts of the instrument are moving.

⚠ CAUTION

Injury due to sharp objects or broken glass

Instrument components, e.g., glass, can break and lead to injuries.

- Always proceed with focus and care.

NOTICE**Damage to the instrument or malfunction due to the use of unsuitable parts**

- Only use parts from METTLER TOLEDO that are intended to be used with your instrument.

NOTICE**Damage to the instrument**

The instrument contains no user-serviceable parts.

- 1 Do not open the instrument.
- 2 In the event of problems, please contact a METTLER TOLEDO representative.

NOTICE**Damage to the instrument due to inappropriate cleaning methods**

If liquid enters the housing, it can damage the instrument. The surface of the instrument can be damaged by certain cleaning agents, solvents, or abrasives.


- 1 Do not spray or pour liquid on the instrument.
- 2 Only use the cleaning agents specified in the Reference Manual (RM) of the instrument or the guide "8 Steps to a Clean Balance".
- 3 Only use a lightly moistened, lint-free cloth or a tissue to clean the instrument.
- 4 Wipe off any spills immediately.

3 Design and Function

3.1 Function description

The **QLX3 liquid module** can be added to **XPR analytical balances** to perform automatic dosing of liquids. The liquid dosing head is attached to the liquid module and located inside the weighing chamber. The vertical position of the liquid module is adjusted manually. The **QL3 pump** and **QLL kit for bottle** are completing the system. The **QL3 pump** is used to build pressure in the bottle. As soon as the pressure is sufficiently high, the micro dispensing valve in the liquid dosing head opens and liquid ascends the liquid tube. Several pumps can be connected to the balance simultaneously and each pump can accommodate up to three bottles (QLL kits).

3.2 Overview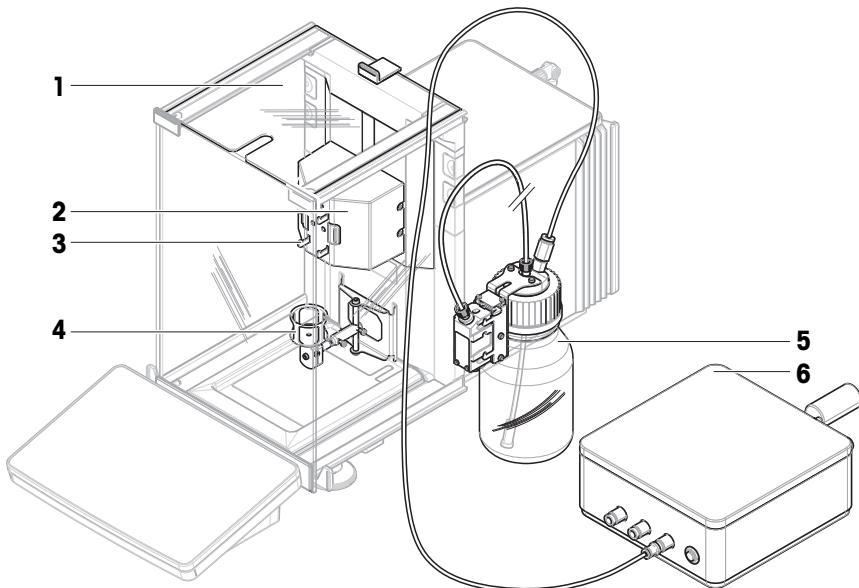


1	Top door	4	ErgoClip vial
2	Liquid module	5	QLL kit for bottle (available separately)
3	Dosing head holder	6	QL3 pump (available separately)

3.3 Components description


Dosing head holder

The dosing head holder is where the liquid dosing head is secured. The dosing head can be released by pressing one of the release buttons on either side of the holder.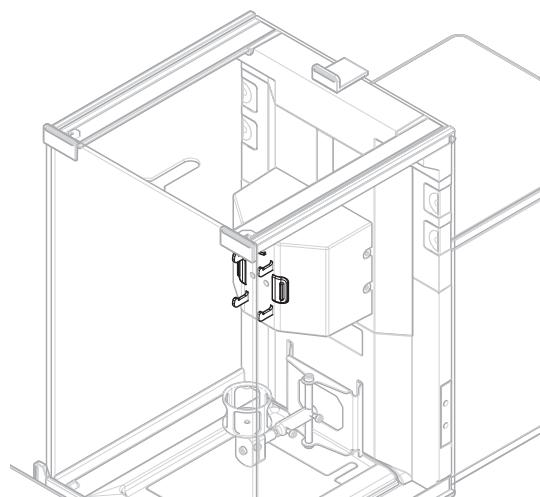


ErgoClip vial

The ErgoClip vial is used together with vial adapters to position vials optimally. Those adapters are available for vials of various diameters and heights. See [Accessories ▶ Page 34].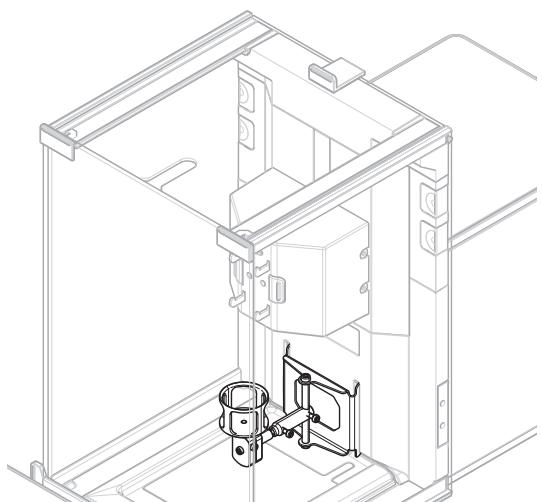


Liquid module

The liquid module is where the liquid dosing head is attached. It can be moved up and down manually to adjust the distance between the sample vessel and the tip of the dosing head.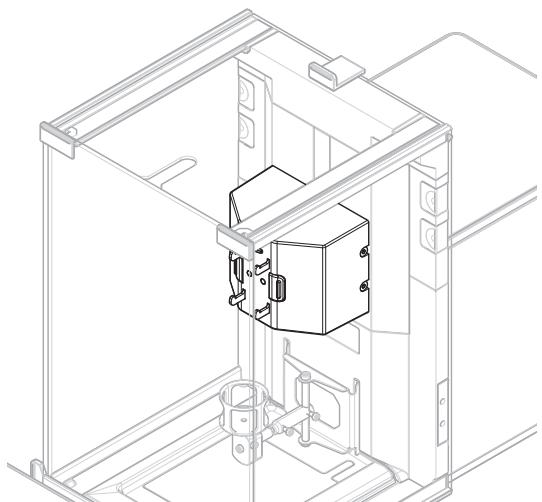

Top door liquid module

The liquid module is provided with a top door designed with a slit through which the liquid tube can pass to reach the inside of the weighing chamber. This door is installed in the same railings as the standard top door of the balance and can be opened and closed manually as well as automatically.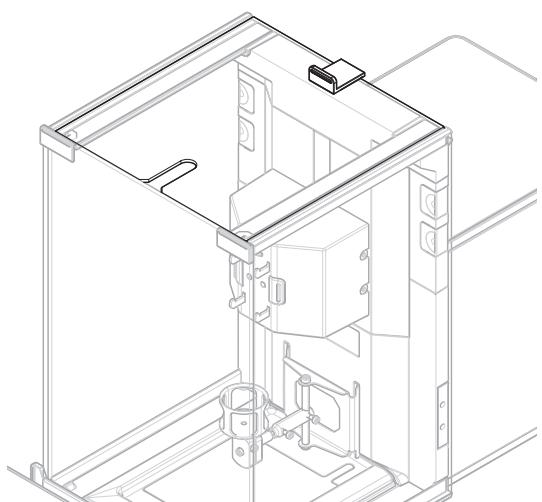


3.4 Dosing head RFID tag

Each powder and liquid dosing head is equipped with an integrated RFID tag (1) that stores and exchanges data with the instrument.

Various data is stored in the RFID tag of the dosing heads, such as the substance name, lot ID, filling date, expiry date, etc. It also includes customized data fields.

This data can be edited on the terminal and should be set before using a new dosing head such that the data is available for reports and labels.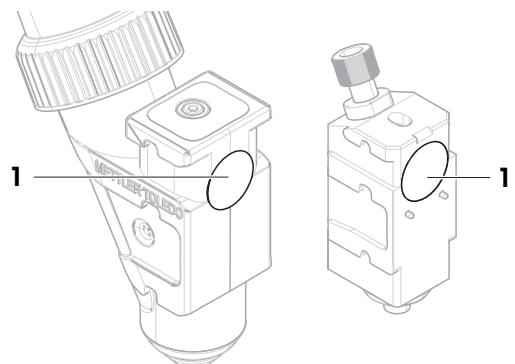

4 Installation and Putting into Operation

This instrument must be installed by a METTLER TOLEDO service technician.

The QLX3 liquid module is compatible with all XPR analytical balances with tall draft shield.

4.1 Scope of delivery

QLX3 liquid module

- QLX3 liquid module
- Top door liquid module
- ErgoClip vial
- Various vial adapters, 4 pcs
- Declaration of Conformity
- User Manual

The scopes of delivery for the **QL3 pump** and the **QLL kits for bottles** are in the documentation delivered with these products.

4.2 Selecting the location


A balance is a sensitive precision instrument. The location where it is placed will have a profound effect on the accuracy of the weighing results.

Requirements of the location


Place indoors on stable table

Avoid direct sunlight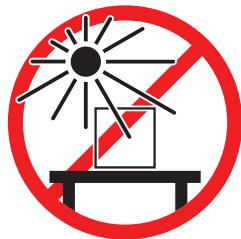


Ensure sufficient spacing

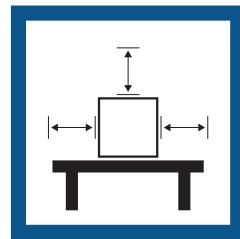
Avoid vibrations


Level the instrument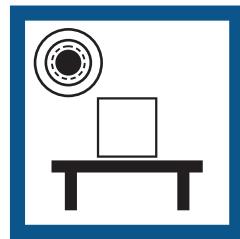

Avoid strong drafts

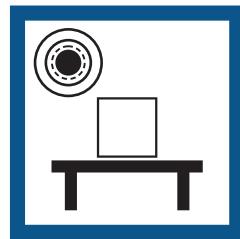Provide adequate lighting

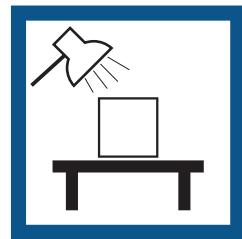Avoid temperature fluctuations


Sufficient spacing for balances: > 15 cm all around the instrument

Take into account the environmental conditions. See "Technical Data".

4.3 Setting up the liquid dosing system

4.3.1 Interface of the pump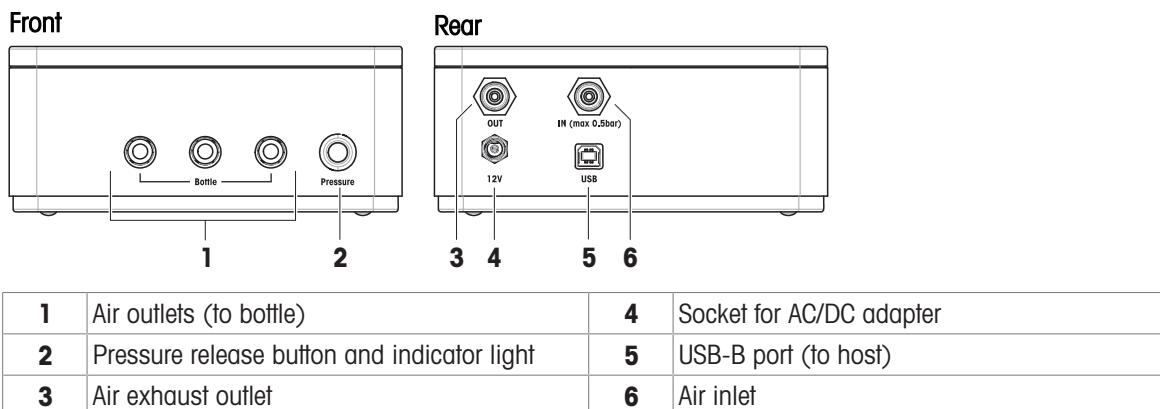

4.3.2 Wiring the pump


- 1 Install the cables in such a way that they cannot be damaged or interfere with operation.
- 2 Insert the plug of the AC/DC adapter in the power inlet of the pump.
- 3 Secure the plug by firmly tightening the knurled nut.
- 4 Insert the plug of the power cable into a grounded power outlet that is easily accessible.
- 5 Using the USB cable, connect the USB-B port of the pump to one of the USB-A ports of the balance.

4.3.3 Connecting the tubes

Tubes definition

The liquid tube is the thinner tube (**1**) used for transporting liquid from the bottle to the liquid dosing head. The air tube is the slightly bigger tube (**2**) used for pumping air into the bottle. By adding air through the air tube, pressure rises in the bottle. When the pressure reaches the target pressure, between 0.3 and 0.5 bar (4.4 to 7.2 psi), the micro dispensing valve in the dosing head opens and liquid can ascend the liquid tube.

- 1** Liquid tube
- 2** Air tube

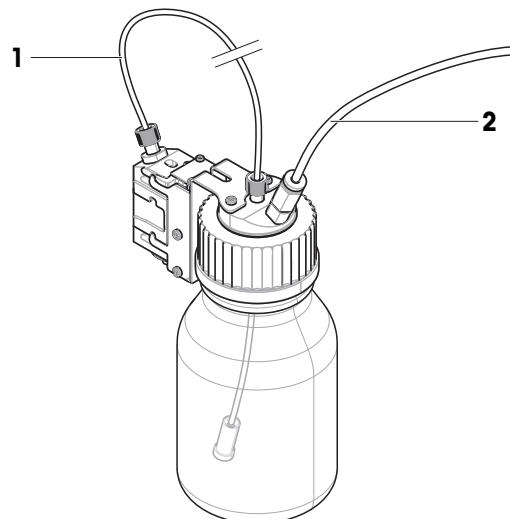Preparing the liquid tube

CAUTION

Injury due to leaking liquids


Wrongly cut tubing can result in leaking connections.

- Cut the tubes with a tube cutter or a sharp knife.

- The liquid dosing head is inserted in the liquid dosing head support of the bottle.

- Using a tube cutter or a sharp knife, cut a sufficient amount of tubing. The appropriate length mainly depends on the distance between the balance and the bottle while dosing.

Recommended length: about 0.9 m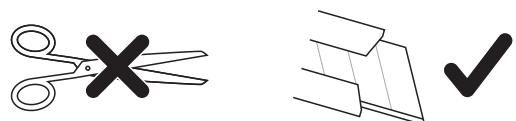

- NOTICE: Leaking liquid due to wrong assembly. Pay attention to the orientation of the sealing ring when threading it on the tube.**

Place the sealing ring (1) on a flat stable surface, e.g., a table or a workbench, with the wider end downwards.

- Take the end of the liquid tube and press it into the sealing ring.


→ This is the dosing-head end of the tube. The opposite end is the bottle end.

- Thread the two fastening nuts (2), paying attention to the orientation.

- NOTICE: Leaking liquid due to wrong assembly. Pay attention to the orientation of the sealing ring when threading it on the tube.**

Thread the sealing ring (3) from the bottle end of the tube.

- Slide the sealing ring until the distance (D) is enough for the tube to reach the bottom of the bottle. Recommended distances for typical bottle volumes are listed next.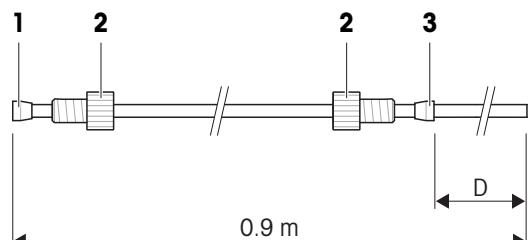


Typical distances (D) between the sealing ring and the bottle-end of the tube

Thread	Bottle volume	Distance (D)
GL45	1000 ml	220 mm
	500 ml	170 mm
	250 ml	135 mm
GL25	25 ml	90 mm

Connecting the liquid tube

- Insert the dosing-head end of the tube (1) in the dosing head (2).
- Tightly fasten the fastening nut to the dosing head.
- Insert the bottle end of the tube through the corresponding hole in the bottle cap (3). The tube should reach the bottom of the bottle.
- If needed, attach the suction filter (4) to the bottle end of the tube.
- Tightly fasten the fastening nut to the bottle cap.
- Screw the cap to the bottle (5).

The suction filter is used to ensure that no particles or impurities are carried through the liquid dosing head. Using the suction filter will extend the lifetime of the dosing head. However, when dosing solutions, molecules of one of the substances might be absorbed by the suction filter, altering the concentration of the solution. The suction filter should only be used when dosing pure solvents.

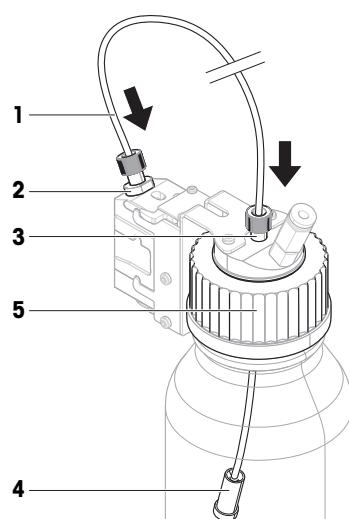Connecting the air tube


WARNING

Injury and/or damage due to reacting substances

When pressure is released from the bottle, the air/gas in the bottle moves back towards the pump. The air/gas coming from the coupled outlets mixes in the pump. Molecules of the substances in the various bottles can get in contact through this contaminated air/gas.

- 1 Do not connect bottles with incompatible liquids to the same pump simultaneously.
- 2 Before connecting a second, incompatible liquid to the pump, disconnect the first bottle and purge the pump cavity with clean air/gas.


- 1 Using a tube cutter or a sharp knife, cut a sufficient amount of tubing. The appropriate length mainly depends on the distance between the bottle and the pump while dosing.
Recommended length: about 0.7 m
- 2 Connect the air tube (1) to the air inlet of the bottle (2)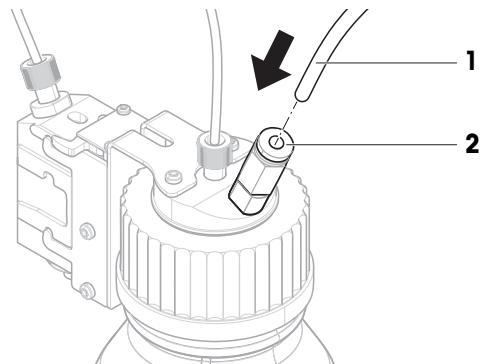

- 3 Insert the other end of the air tube (1) in a tube adapter (3) and press firmly.
- 4 Connect the tube adapter (3) to one of the air outlets (4) of the pump. Press until you hear a click.


When a tube is connected to the air outlet of the pump, the valve of the air outlet opens. Never leave a tube that is connected to the air outlet unconnected at the other end because pressure cannot build up.

You can connect up to three bottles to each pump.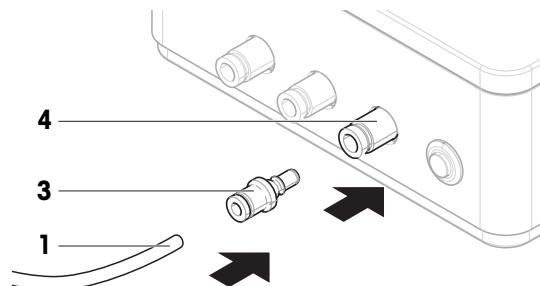

Connecting the muffler

- Insert the muffler (1) into the air inlet (2) to absorb the noise.

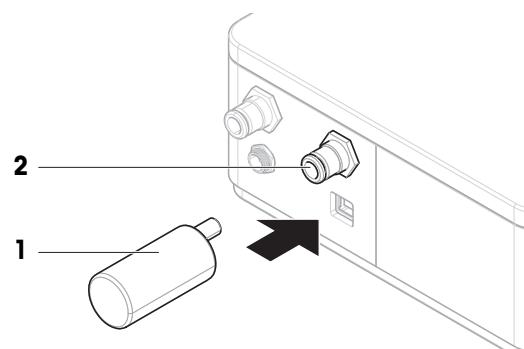Removing the air tube

NOTICE


Damage to tube connectors due to mishandling

If the tubes are not removed correctly, the connectors of the pump and the bottle cap can be damaged.



- To remove the tubes, press down the ring on the connector and pull out the tube carefully.

- 1 Remove the tube adapter (1) from the pump by pushing the connector (2) of the air outlet towards the pump.

→ The adapter is released and can be removed.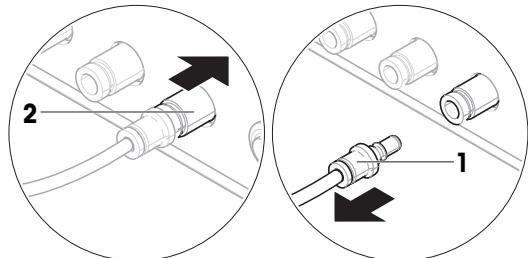

- 2 Remove the air tube from the bottle by firmly pressing down the ring (3) and pulling the tube (4) simultaneously.
- 3 If necessary, e.g., for maintenance purposes, use the same procedure to remove the air tube from the tube adapter: firmly press the ring and pull the tube simultaneously.

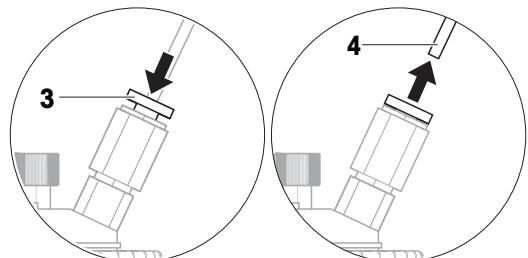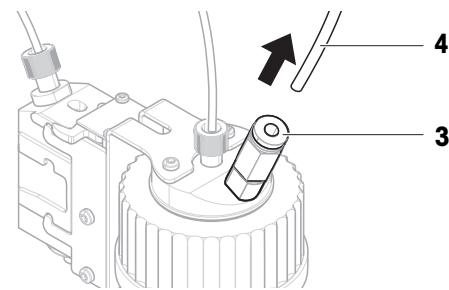Using the pump with external gas

The liquid can be protected by feeding an external gas, e.g., nitrogen, to the pump. Make sure that the pressure of the external gas does not exceed 0.5 bar (7.2 psi).

WARNING

Injury and damage to the pump or the bottle due to high pressure

High pressure from external gas can damage the pump or the bottle.


- 1 Use a regulator on the external gas line.
- 2 Ensure that the pressure of the external gas does not exceed 0.5 bar (7.2 psi).

- A regulator is connected to the external gas line.
- 1 Remove the muffler from the air inlet (1).
 - 2 Connect the tube of external gas (2) to the air inlet (1).

Note

Outer tube diameter: 6 mm

Pressure in external gas line: The pressure must be at least 0.1 bar (1.5 psi). The pressure must not exceed the configured dosing pressure set on the terminal. The external gas tube is not provided by METTLER TOLEDO.

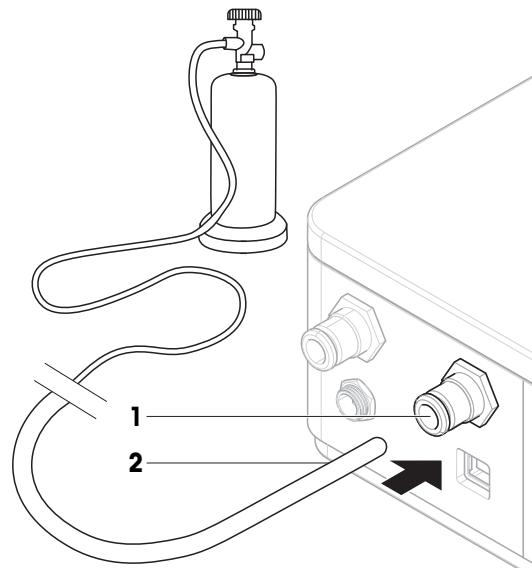Collecting contaminated air

WARNING

Injury or death due to toxic, explosive, or flammable substances

If you use toxic, explosive, or flammable liquids with the pump, the exhaust air will be contaminated.


- Connect a tube to the exhaust air outlet to collect the contaminated air.

- Connect a tube (1) to the exhaust air outlet (2) to collect the contaminated air into a safe container (3).

Note

Outer tube diameter: 6 mm

The exhaust air tube and the container are not provided by METTLER TOLEDO.

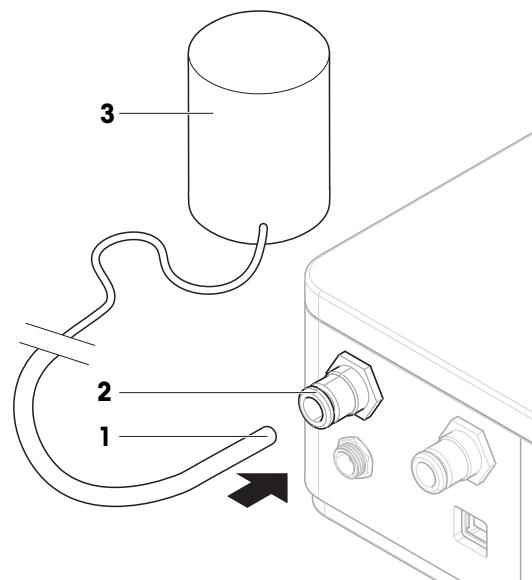5 Operation

5.1 Liquid dosing


For further information, consult the Reference Manual (RM) of your XPR balance.

► www.mt.com/XPR-analytical-RM

5.1.1 Installing and removing the liquid dosing head


Installing the dosing head

- 1 Slide the dosing head (1) onto the dosing head holder (2) until it comes to a stop.
- 2 Slightly press down the dosing head until it properly sits in the dosing head holder (2).
 - ⇒ The dosing head is locked in the holder.
- 3 Thread the liquid tube through the slot in the top door.
 - ⇒ The dosing head is ready to dose.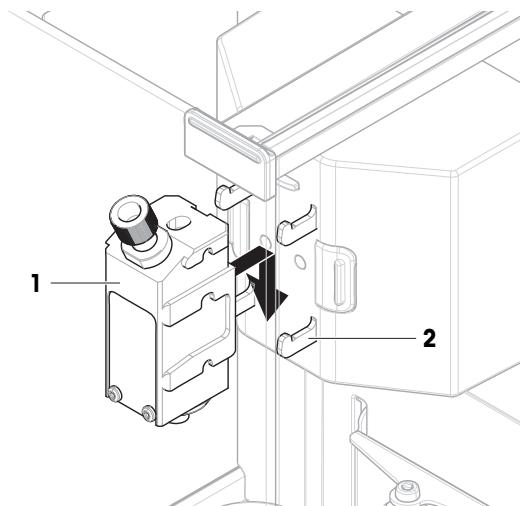

Removing the dosing head

- 1 Press the release button (1) on either side of the dosing head holder to release the dosing head (2).
- 2 Simultaneously, pull the dosing head (2) up and outwards.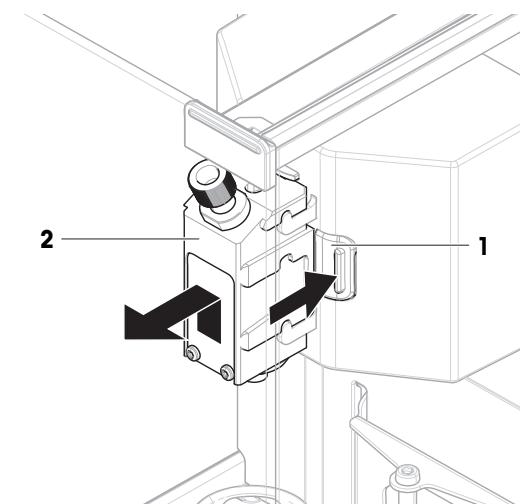


5.1.2 Adjusting the dosing height


The vertical position of the liquid module is adjusted by moving it manually up and down along its rails.

5.1.3 Attaching the dosing head to the bottle cap

- 1 Insert the liquid dosing head in the liquid dosing head support.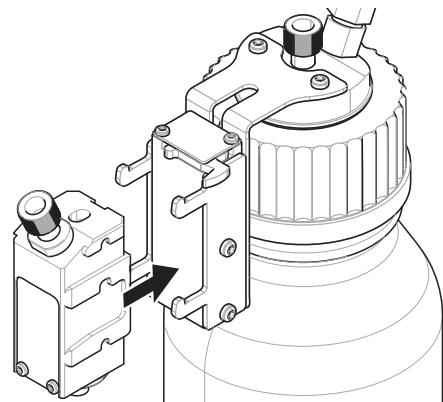


- 2 To remove the liquid dosing head from the liquid dosing head support, pull the lever (1) towards the dosing head and remove the liquid dosing head (2).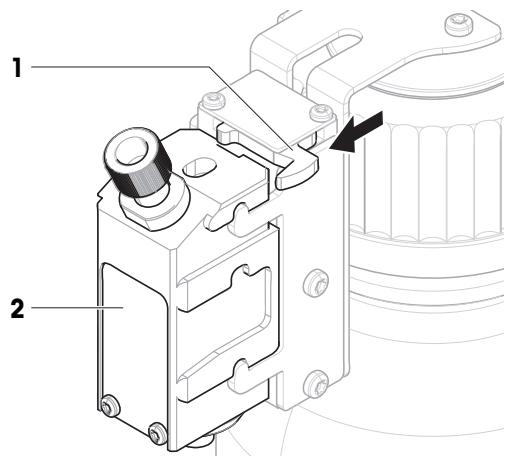


5.1.4 Handling the bottle

Filling the bottle

The suction filter (1) has to be covered with liquid at all times. Before the suction filter gets dry, refill the bottle.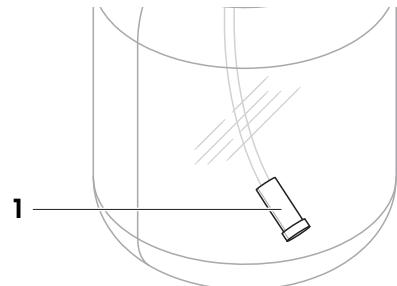

- The pressure is released.
- 1 **⚠ CAUTION: Injury due to splashing liquids. Make sure that the pressure in the bottle is released.**
Unscrew the cap.
 - 2 Fill in the liquid. Do not exceed the maximum (max. is shown on bottle, e.g., 1000 ml). The air above the liquid is necessary for dosing.
 - 3 Screw the cap on tightly.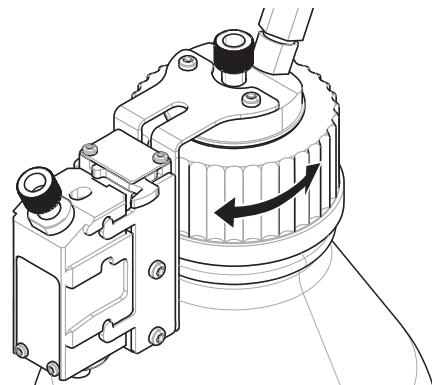

Changing the bottle content

This procedure is used if you have a single cap with dosing head and want to change the liquid to be dosed.

- The pressure is released.


- 1 **⚠ CAUTION: Injury due to splashing liquids. Make sure that the pressure in the bottle is released.**
Unscrew the cap.
- 2 Remove the suction filter, if applicable.
- 3 If the cap needs to be cleaned, remove the air tube.
Unscrew the fastening nut from the cap.
Rinse the cap with the appropriate solvent or liquid.
Insert the liquid tube into the cap.
- 4 If the liquid tube needs to be cleaned with a solvent, fill the bottle with the appropriate solvent.
Screw the cap on the bottle.
Insert the air tube on the cap.
Purge using the **Purge** function on the terminal.
Unscrew the cap.
Dispose of the remainder of solvent.
- 5 Attach a new suction filter, if applicable.
- 6 Screw the cap on the bottle containing the new liquid.
- 7 Check that the cap is tight.
- 8 Connect the air tube to the new bottle.
- 9 Purge using the **Purge** function.

Connecting the air tube to a different bottle

If you have more than one bottle equipped with cap and dosing head, and you want to use the same pump outlet and air tube to dose from another bottle:

- The pressure is released.

- 1 Install the dosing head on the dosing head support of the bottle.
- 2 Unplug the air tube by pressing down the ring (1) and simultaneously pulling out the tube (2).
- 3 To seal the bottle, insert the pin delivered with the QLL kit (3) into the air tube connector.
- 4 Take the new bottle.
- 5 Connect the air tube to the new bottle.
- 6 To continue dosing with the new bottle, install the dosing head.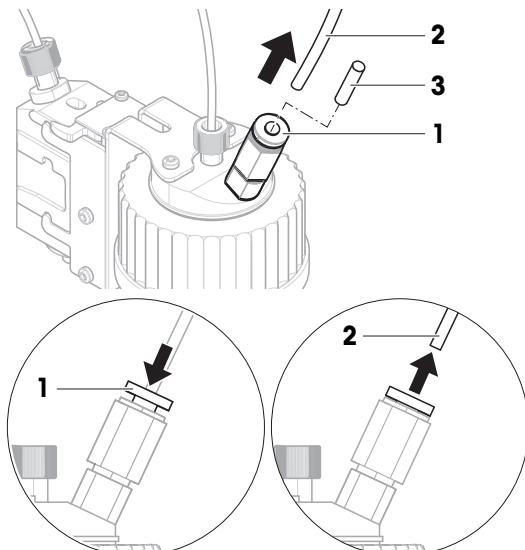

5.1.5 Using the QL3 pump


Pressure indicator light

The pressure indicator light displays the status of the pump:

- Light on: pressure is building up or pressure is established
- Light off: there is no pressure and no pressure is building up
- Light blinking: pump error and/or warning

Releasing the pressure

- The pressure indicator light (**1**) is on.
- Press the pressure release button (**2**) to release the pressure.
- The status light (**1**) turns off when pressure is released.

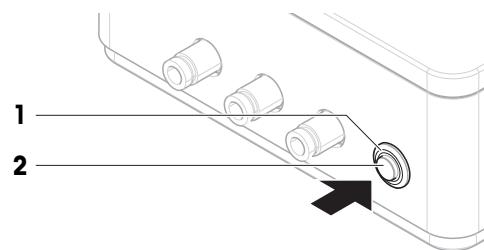Purging the pump cavity

If bottles containing incompatible liquids (and for which the fumes should not mix) are connected to the pump one after the other, it is recommended to purge the pump cavity before connecting the second bottle to the pump.

- A dosing head is attached to the dosing or liquid module. The bottle of this dosing head is not connected to any pump.
 - A method **Automated dosing** or **Automated solution prep.** is running on the terminal.
- 1 Disconnect all tube adapters from the pump.
 - 2 Connect an empty tube adapter to the right-most air outlet at the front of the pump.
 - The pump is trying to build pressure and air flows through the cavity of the pump, purging it.
 - The pump cavity has been purged and bottles can safely be re-connected to the air outlets.

5.2 Using other weighing methods

To use other methods with the balance, move the liquid module to its uppermost position.

For further information, consult the Reference Manual (RM) of your XPR balance.

► www.mt.com/XPR-analytical-RM

6 Maintenance

The appropriate maintenance interval depends on your standard operating procedure (SOP).

Please contact your METTLER TOLEDO representative for details about the available service options. Regular servicing by an authorized service technician ensures constant accuracy for years to come and prolongs the service life of your instrument.

6.1 Cleaning

NOTICE

Damage to the instrument due to inappropriate cleaning methods

If liquid enters the housing, it can damage the instrument. The surface of the instrument can be damaged by certain cleaning agents, solvents, or abrasives.

- 1 Do not spray or pour liquid on the instrument.
- 2 Only use the cleaning agents specified in the Reference Manual (RM) of the instrument or the guide "8 Steps to a Clean Balance".
- 3 Only use a lightly moistened, lint-free cloth or a tissue to clean the instrument.
- 4 Wipe off any spills immediately.

For further information on cleaning a balance, consult "8 Steps to a Clean Balance".

► www.mt.com/lab-cleaning-guide

Detailed information about the compatibility of cleaning agents can be found in the Reference Manual (RM) of your XPR balance.

► www.mt.com/XPR-analytical-RM

6.1.1 Cleaning the housing

The housing material of the liquid module is the same as the one of the balance. All surfaces can therefore be cleaned with a commercially available, mild cleaning agent.

6.1.2 Cleaning the weighing chamber

The cleaning procedure is the same as for the balance. Refer to the Reference Manual of the balance for more details.

6.1.3 Cleaning the liquid dosing elements

Cleaning the suction filter

- 1 **Once a week**, check visually that the suction filter is clean. If necessary, rinse the suction filter with solvent using the **Purge** function on the terminal or replace the suction filter.
- 2 At least **once a year**, change suction filter. Maintenance interval depends on the liquid used.

Purging the liquid dosing head

The liquid dosing head can be rinsed/purged by letting a large quantity of solvent (or another liquid) run through it. Use the **Purge** function. See the Reference Manual of the XPR balance.

- The bottle is filled with enough solvent to purge the dosing head.
- A big sample vessel is on the weighing pan to collect the solvent used to purge the dosing head.
- The liquid dosing head is installed on the dosing or liquid module.
- Use the **Purge** function on the terminal to rinse the dosing head.


Cleaning the micro dosing valve

The liquid dosing heads QL001 have a micro dosing valve that can be disassembled and cleaned, for example, in an ultrasonic bath.

Note

The QL003 dosing head does not contain a micro dosing valve. Cleaning this dosing head using the **Purge** function is sufficient (see above).


- The liquid dosing head is installed on the dosing or liquid module.
 - A sample vessel is on the weighing pan, big enough for the quantity of liquid in the liquid tube.
 - The pressure is released.
- 1 **⚠ CAUTION: Injury due to splashing liquids. Make sure that the pressure in the bottle is released.**
Empty the liquid tube by replacing the bottle by an empty bottle and using the **Purge** function.
→ The liquid tube is empty.
 - 2 Release pressure.
 - 3 **⚠ CAUTION: Injury due to splashing liquids. Make sure that the pressure in the bottle is released.**
Remove the dosing head from the dosing or liquid module.
 - 4 **⚠ WARNING: Injury or contamination from hazardous substances. Be aware of any liquid draining from the tubing and the valve.**
Open the dosing head with the micro dosing valve tool (1), if applicable.
 - 5 Take out the micro dosing valve and clean it, e.g., in an ultrasonic bath.
 - 6 After the cleaning, reinstall the micro dosing valve and refill/replace the bottle.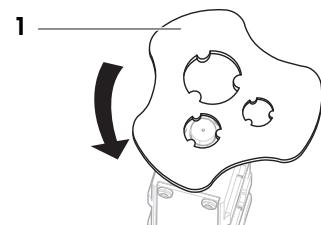


6.2 Replacing the sealing ring and fastening nut on the liquid tube

Removing the sealing ring and the fastening nut

If liquid is leaking from the dosing head nut, replace the sealing ring and fastening nut at the dosing head (1). If pressure cannot be built in the bottle, replace the sealing ring and fastening nut at the bottle cap (2).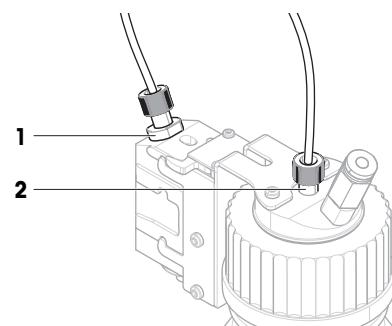

- The pressure is released.
- 1 **⚠️ CAUTION: Injury due to splashing liquids. Make sure that the pressure in the bottle is released.**
Unscrew the fastening nut at the dosing head or the bottle cap.
 - 2 To access the sealing ring (1), slide back the fastening nut (2).
 - 3 If changing the nut and ring at the bottle cap, remove the suction filter and slide the sealing ring all the way to the bottle-end of the tube.
 - 4 Using a tube cutter or a sharp knife, cut off the tube above the sealing ring (1).
 - 5 Remove the fastening nut.

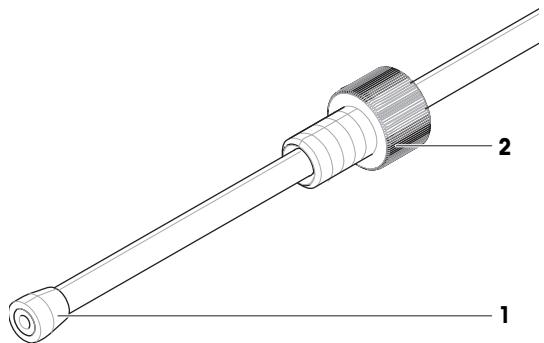Reconnecting the liquid tube

Insert the new fastening nut sealing ring, and reconnect the liquid tube.

See also

📄 Connecting the tubes ▶ Page 14

7 Troubleshooting

For further information, consult the Reference Manual (RM) of your XPR balance.

► www.mt.com/XPR-analytical-RM

7.1 Error symptoms

Error symptom	Possible cause	Diagnostic	Remedy
The liquid dosing head does not dispense any liquid.	The suction filter is clogged. The dosing head is not installed properly.	Check if there is liquid in the liquid tube. –	Clean or replace the suction filter. Remove and reinstall the dosing head in the holder. Make sure to press until you hear a click.
Liquid leaks from the bottle cap and/or the dosing head.	The tube is not properly connected.	–	Tighten the fastening nut at the bottle cap and/or the dosing head. Make sure that the end of the tube is cut straight.
	The sealing ring and/or fastening nut are damaged.	–	Change the sealing ring and the fastening nut at the bottle cap and/or the dosing head. See "Maintenance". Make sure that the end of the tube is cut straight.
Liquid drips from the QL003 liquid dosing head.	Impurities have entered the QL003 liquid dosing head.	–	Use the Purge function for at least 10 seconds to rinse the dosing head. Check if the dosing head is still dripping. Repeat if necessary. If the dosing head cannot be cleaned properly, replace it.
The status light of the pump does not blink when the device is connected to power.	The pump is disconnected from power.	Disconnect and reconnect the AC/DC adapter from the pump. The indicator light should blink once when the device is connected to power. Check that the AC/DC adapter and the power cable are not damaged.	Replace the AC/DC adapter and power cable.
	The pump is damaged.	Check with another pump, if available.	Replace the pump. Contact your METTLER TOLEDO service representative.

Error symptom	Possible cause	Diagnostic	Remedy
On the display, the QL3 pump does not appear in the list of devices connected to the balance.	The pump is disconnected from power.	Disconnect and reconnect the AC/DC adapter from the pump. The indicator light should blink once when the device is connected to power. Check that the AC/DC adapter and the power cable are not damaged.	Replace the AC/DC adapter and power cable.
	The USB cable is not properly connected.	Check that the USB cable is connected properly.	Connect the USB cable properly.
	The USB cable is damaged.	Check that the USB cable is not damaged.	Replace the USB cable.
	The USB-A port of the balance is damaged.	Disconnect the pump from the USB-A port of the balance. Connect a USB mouse to the same USB-A port. Verify that a pointer (arrow) appears on the terminal and can be moved by moving the mouse.	If the mouse pointer does not appear, contact your METTLER TOLEDO service representative.
	The pump is damaged.	Check with another pump, if available.	Replace the pump. Contact your METTLER TOLEDO service representative.

8 Technical Data

8.1 General data

Weight (without packaging):	470 g
Power consumption:	12 V DC ± 6%, 1 A

Protection and standards

Overvoltage category:	II
Degree of pollution:	2
Range of application:	Use only indoors in dry locations

Environmental conditions

Height above mean sea level:	Up to 5000 m
Ambient temperature:	+5 – +40 °C
Relative air humidity:	20% to max. 80% at 31 °C, decreasing linearly to 50% at 40 °C, non-condensing

Storage conditions (in packaging)

Ambient temperature:	-25 – +70 °C
Relative air humidity:	10 – 90%, non-condensing

8.2 Model-specific data

Liquid dosing heads

	QL001	QL003
Dosing properties		
Suitable liquids: max. viscosity ¹⁾	20 mPa·s	20 mPa·s
Typical values²⁾		
Repeatability (sd)	1 mg	2 mg
Dosing time ³⁾	40 s	25 s

¹⁾ 1 mPa·s = 1 cP = 1 cSt = 1 mm²/s

²⁾ for H₂O

³⁾ for 10 g

Tubing

	Outer diameter	Inner diameter	Recommended length
Liquid tube, for GL45 bottles (dosing head to bottle)	3.2 mm	1.6 mm	0.9 m
Liquid tube, for GL25 bottles (dosing head to bottle)	1.6 mm	0.8 mm	0.9 m
Air tube (pump to bottle)	4.0 mm	2.4 mm	0.7 m
Tube for exhaust air	6 mm	–	–
Tube for external gas	6 mm	–	–

Suction filter

Tubing outer diameter	3.2 mm
Filter pore size	10 µm

Bottle

Pressure resistance, min. ¹⁾	1.5 bar
Burst pressure, min.	3 bar
Volume, max.	2 l

¹⁾ According to DIN EN 1595: Pressure Equipment made from Borosilicate Glass 3.3 – General Rules for Design, Manufacture and Testing

QL3 pump

Maximum pressure ¹⁾	1.5 bar
--------------------------------	---------

¹⁾ In the unlikely event of a malfunction, the overpressure valve opens if the maximum pressure is reached, to preserve the integrity of the system.

8.3 Material specifications

Materials in contact with the substance to be dosed. Please note that traces of all contact material could be transferred to the sample.

8.3.1 Glossary of material acronyms

Acronyms defined in ISO 1043: Plastics – Symbols and abbreviated terms.

ETFE	= Ethylene tetrafluoroethylene
FEP	= Fluorinated ethylene propylene
FFKM	= Perfluoroelastomer
PE	= Polyethylene
PEEK	= Polyetheretherketone
PMP	= Polymethylpentene
POM	= Polyoxymethylene
PP	= Polypropylene
PP (ESD)	= Polypropylene (electrostatic dissipative)
PTFE	= Polytetrafluoroethylene

8.3.2 Liquid dosing head QL001

	Designation	Material
1	Fastening nut	PEEK
2	Sealing ring	ETFE
3	Tube connector	Stainless steel 1.4404
4	O-Ring	FFKM
5	Solenoid valve	PEEK Sapphire-Ruby Stainless steel 1.4105IL Stainless steel 1.14301 / 1.4306 Stainless steel 1.4305

8.3.3 Liquid dosing head QL003

	Designation	Material
1	Fastening nut	PEEK
2	Sealing ring	ETFE
3	Nozzle	Ceramic Al ₂ O ₃ > 99.7%
4	Valve block	PTFE
5	Gasket	EPDM
6	Valve	PEEK

8.3.4 QLL standard/advanced kit for bottles


The QLL kits for bottle include a liquid dosing head, see [Liquid dosing head QL001 ▶ Page 30], [Liquid dosing head QL003 ▶ Page 31].

	Designation	Material
1	Cap insert	PE
2	Bottle	Borosilicate glass
3	Tubing	FEP
4	Suction filter	PP

8.4 Dimensions

QLX3 liquid module and XPR balance

Dimensions in mm.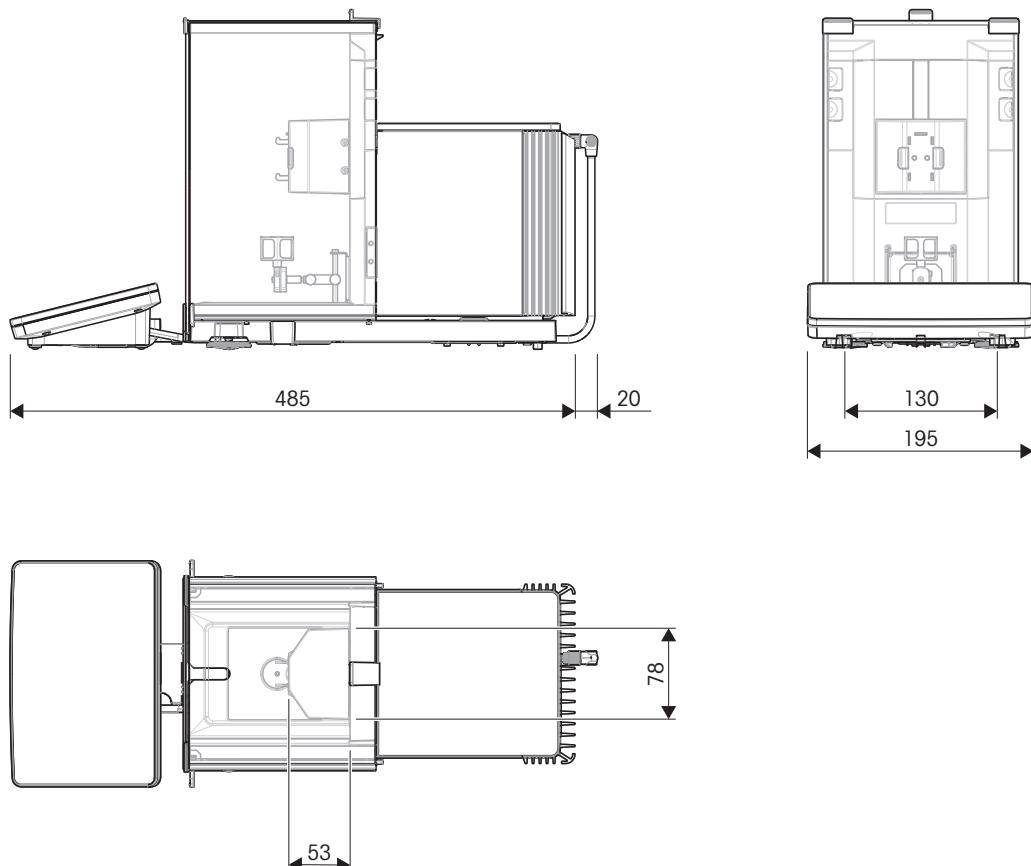


Dimensions

Maximum height of sample vessel, without adapter, when placed on SmartGrid weighing pan	140 mm
Minimum opening of sample vessel (diameter)	6 mm
Vertical range of movement of liquid module	53 mm

QLX3 liquid module

Dimensions in mm.

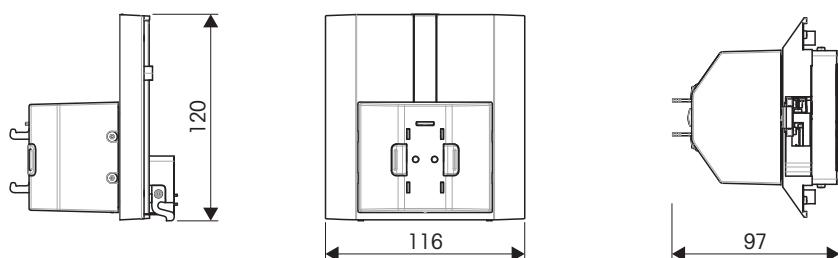9 Disposal

In conformance with the European Directive 2012/19/EU on Waste Electrical and Electronic Equipment (WEEE) this device may not be disposed of in domestic waste. This also applies to countries outside the EU, per their specific requirements.

Please dispose of this product in accordance with local regulations at the collecting point specified for electrical and electronic equipment. If you have any questions, please contact the responsible authority or the distributor from which you purchased this device. Should this device be passed on to other parties, the content of this regulation must also be related.

10 Accessories and Spare Parts

10.1 Accessories

Accessories are additional components that could help you in your workflow.

Please contact your METTLER TOLEDO sales representative for more details about using accessories with your product.

	Description	Order no.
Printers		
	CLS-631 label printer (RS232C/USB-A) Label and ink ribbon kit The NetCom Kit might be needed for this printer. Contact your local METTLER TOLEDO service representative.	11141820 30004309
	P-52RUE dot matrix printer RS232C, USB and Ethernet connections, simple print-outs Paper roll (length: 20 m), set of 5 pcs Paper roll (length: 13 m), self-adhesive, set of 3 pcs Ribbon cartridge, black, set of 2 pcs	30237290 00072456 11600388 00065975
	P-56RUE thermal printer with RS232C, USB and Ethernet connections, simple print-outs, date and time Paper roll, white (length: 27 m), set of 10 pcs Paper roll, white, self-adhesive (length: 13 m), set of 10 pcs	30094673 30094723 30094724
	P-58RUE thermal printer with RS232C, USB and Ethernet connections, simple print-outs, date and time, label printing, balance applications, e.g., statistics, formulation, totaling Paper roll, white (length: 27 m), set of 10 pcs Paper roll, white, self-adhesive (length: 13 m), set of 10 pcs Paper roll, white, self-adhesive labels (550 labels), set of 6 pcs Dimension of the label 56×18 mm	30094674 30094723 30094724 30094725
ErgoClips		
	ErgoClip adapter holder	30521809
	ErgoClip height extension	30542824

	ErgoClip adapter stabilizer	30542831
	ErgoClip vial	30521808

RFID readers / writers / cards

	EasyScan USB Reads and writes RFID tags.	30416173
	Smart Tag Set of 50 pieces Set of 200 pieces	30101517 30101518

Barcode readers

	Corded USB barcode reader	30417466
---	---------------------------	----------

Adapters

	Vial adapters (POM), diameter < 15 cm 8.5 mm × 15 mm (5 pcs.)	30428901
	9 mm × 25 mm (5 pcs.)	30428902
	9.5 mm × 25 mm (5 pcs.)	30428903
	10 mm × 25 mm (5 pcs.)	30428904
	10.5 mm × 25 mm (5 pcs.)	30428905
	11 mm × 30 mm (5 pcs.)	30428906
	11.5 mm × 30 mm (5 pcs.)	30428907
	12 mm × 20 mm (5 pcs.)	30428908
	12.5 mm × 30 mm (5 pcs.)	30428909
	13.5 mm × 20 mm (5 pcs.)	30428910
	14.5 mm × 25 mm (5 pcs.)	30428911

**Vial adapters (POM), diameter > 15 cm**

15.5 mm × 25 mm (5 pcs.)	30428912
16.5 mm × 25 mm (5 pcs.)	30428913
17.5 mm × 25 mm (5 pcs.)	30428914
18.5 mm × 25 mm (5 pcs.)	30428915
19.5 mm × 25 mm (5 pcs.)	30428916
20.5 mm × 25 mm (5 pcs.)	30428917
21.5 mm × 25 mm (5 pcs.)	30428918
22.5 mm × 25 mm (5 pcs.)	30428919
23.5 mm × 25 mm (5 pcs.)	30428920
26 mm × 25 mm (5 pcs.)	30428926
27.5 mm × 25 mm (5 pcs.)	30428921
28.5 mm × 25 mm (5 pcs.)	30428922
29.5 mm × 35 mm (5 pcs.)	30428923
31.5 mm × 30 mm (5 pcs.)	30428924

**Bottle adapters (POM)**

33.5 mm × 15 mm (5 pcs.)	30459921
35.5 mm × 15 mm (5 pcs.)	30459922
37.5 mm × 15 mm (5 pcs.)	30459923
40.5 mm × 15 mm (5 pcs.)	30459924
44.5 mm × 15 mm (5 pcs.)	30459925
48 mm × 15 mm (5 pcs.)	30459926
52 mm × 15 mm (5 pcs.)	30459927

**Tube adapters (POM)**

Eppendorf 1.5 ml (5 pcs.)	30306209
Falcon 16.4 mm × 60 mm (5 pcs.)	30459992
Falcon 29.3 mm × 60 mm (5 pcs.)	30459991

Dosing components

QL3 pump

30418660

QLL standard kit for bottles


- QL001 liquid dosing head
- 250 ml bottle, pressure resistant
- Bottle cap (GL45) with tube connector and dosing head support
- Spare parts for QLL standard kit

30008318

	QLL standard kit for small bottles <ul style="list-style-type: none"> • QL001 liquid dosing head • 25 ml bottle, pressure resistant • Stand for bottle • Bottle cap (GL25) with tube connector • Spare parts for QLL standard kit small 	30237340
	QLL advanced kit for bottles <ul style="list-style-type: none"> • QL003 liquid dosing head • 250 ml bottle, pressure resistant • Bottle cap (GL45) with tube connector and dosing head support • Spare parts for QLL standard kit 	30521817
	Holder for 250 ml bottles	30542827
Various		
	Cable box	11141845
	EasyHub USB	30468768
	Drip pan, gray	30460856


10.2 Spare parts

10.2.1 QLX3 liquid module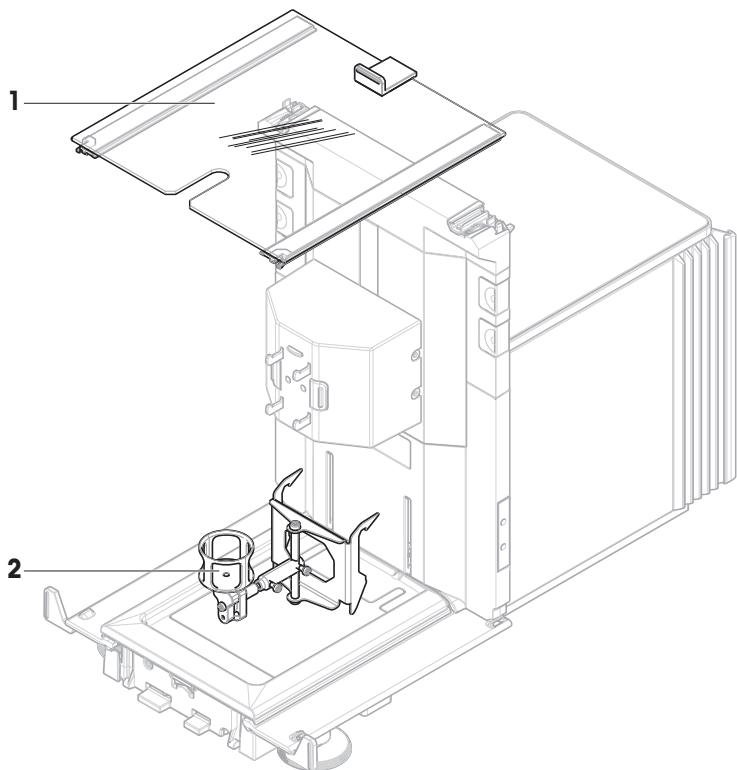


	Order no.	Designation	Remarks
1	30525849	Door top draft shield liquid module	Material: glass
2	30521808	ErgoClip vial	–

10.2.2 QL3 pump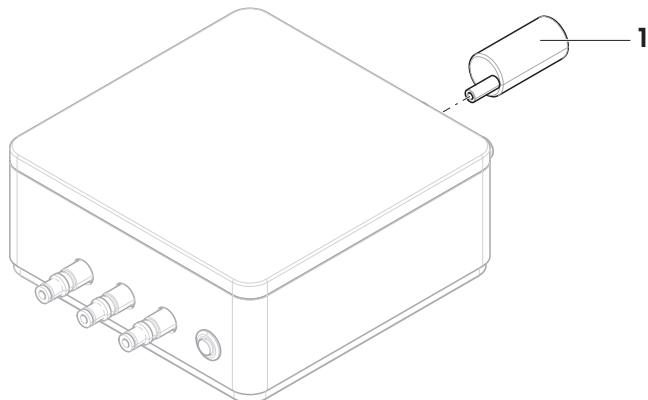

	Order no.	Designation	Remarks
1	30378287	Muffler for pump	Including: muffler, adapter

10.2.3 QLL spare parts kit for bottles

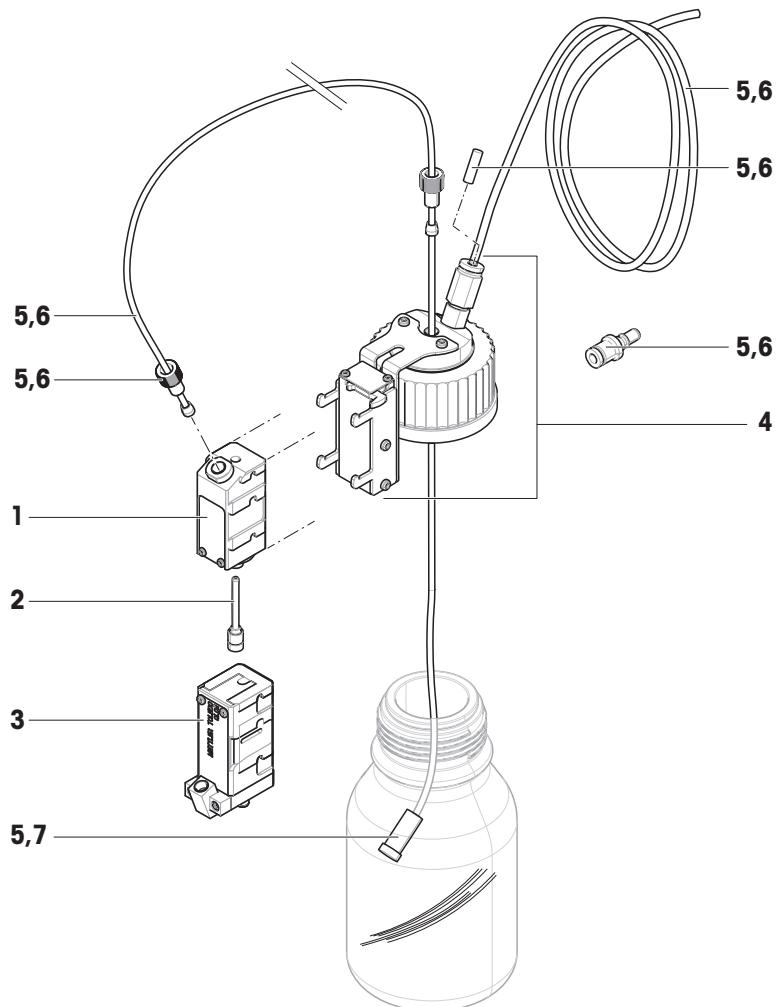	Order no.	Designation	Remarks
1	30080473	Dosing head liquid QL001	–
2	30348196	Micro dosing valve	Including: micro dosing valve tool
3	30542814	Dosing head liquid QL003	–
4	30080708	Bottle cap with tube connector and support	Thread: GL45
5	30025649	Spare parts for QLL kit	Including: tubing OD 4 mm, ID 2.4 mm (2 m), tubing OD 3.2 mm, ID 1.6 mm (2.5 m), 10 sealing rings for tubing OD 3.2 mm, 5 fastening nuts for tubing OD 3.2 mm, 2 closure pins ISO 2338 4H8x20 A1, 2 suction filters for tubing OD 3.2 mm, tube adapter
6	30460218	Spare parts for QLL kit small	Including: tubing OD 4 mm, ID 2.4 mm (2 m), tubing OD 1.6 mm, ID 0.8 mm (2.5 m), 10 sealing rings for tubing OD 1.6 mm, 5 fastening nuts for tubing OD 1.6 mm, 2 closure pins ISO 2338 4H8x20 A1, tube adapter
7	30542832	Suction filter for tubing OD 3.2 mm	Including: 2 suction filters

Index

A

altitude 28

B

bottle

- change 21
- change content 21
- dosing head 20
- fill 20

C

change

- bottle 21
- liquid 21

clean

- housing 23
- liquid dosing head 24
- suction filter 23

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© Mettler-Toledo GmbH 01/2022