InPro 5000 (i) CO₂ Sensors

**Instruction manual
Bedienungsanleitung
Instructions d'utilisation**

InPro 5000 (i)
52 002 465 F

English **Page 3**

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InPro 5000 (i) CO₂ Sensors

Instruction manual

InPro, InFit, ISM and InTrac are trademarks
of the METTLER TOLEDO Group.

Contents

1	Introduction	5
2	Important notes	6
2.1	Notes on operating instructions	6
2.2	Intended use	6
2.3	Safety instructions.....	7
2.4	Environmental protection	7
3	Product description	8
3.1	General information.....	8
3.2	Principles of CO ₂ measurement	8
3.3	Scope of delivery	8
3.4	Equipment features	9
4	Preparation	10
4.1	Start-up	10
4.2	Calibration	11
4.2.1	Connecting the InPro 5000 (i) to cable	11
4.2.2	Connecting the cable to the transmitter	12
4.2.3	Calibration of pH electrode	12
4.3	Mounting the membrane body.....	13
4.4	Mounting the sensor.....	14
4.5	Sterilization	14
5	Operation	15
5.1	Calibration of the CO ₂ sensor.....	15
6	Maintenance	16
6.1	Inspection	16
6.1.1	Visual inspection	16
6.1.2	Checking wiring and transmitter with the pH simulator	16
6.2	Changing the CO ₂ electrolyte and membrane body.....	16
6.3	Maintenance of the interior body (pH electrode)....	17
6.4	Replacement of the interior body (pH electrode) ...	17
6.5	ISM	17
7	Storage	18
8	Product specification	18
8.1	Certificates	18
8.2	Specifications	18
9	Ordering information	19
9.1	Sensors	19
9.2	Accessories.....	19
9.3	Spare parts	19
9.4	Transmitter	20

1 Introduction

Thank you for buying the **InPro™ 5000 (i) sensor from METTLER TOLEDO.**

The construction of the InPro 5000 (i) sensor employs leading edge technology and complies with safety regulations currently in force. Notwithstanding this, improper use could lead to hazards for the user or a third-party, and / or adverse effects on the plant or other equipment.

Therefore, the operating instructions must be read and understood by the persons involved before work is started with the sensor.

The instruction manual must always be stored close at hand, in a place accessible to all people working with the InPro 5000 (i).

If you have questions, which are not or insufficiently answered in this instruction manual, please contact your METTLER TOLEDO supplier. They will be glad to assist you.

2 Important notes

2.1 Notes on operating instructions

These operating instructions contain all the information needed for safe and proper use of the InPro 5000 (i) sensor.

The operating instructions are intended for personnel entrusted with the operation and maintenance of the sensors. It is assumed that these persons are familiar with the equipment in which the sensor is installed.

Warning notices and symbols

This instruction manual identifies safety instructions and additional information by means of the following symbols:

This symbol draws attention to **safety instructions and warnings of potential danger** which, if neglected, could result in injury to persons and / or damage to property.

This symbol identifies **additional information and instructions** which, if neglected, could lead to defects, inefficient operation and possible loss of production.

2.2 Intended use

InPro 5000 (i) sensors are intended solely for in line measurement of dissolved CO₂ (carbon dioxide), as described in this instruction manual.

Any other use, or use not mentioned here, that is incompatible with the technical specifications is deemed inappropriate. The operator is solely responsible for any damage arising from such use.

Other prerequisites for appropriate use include:

- observing the instructions, notes and requirements set out in this instruction manual.
- observing all local safety regulations concerning safety on work.
- observing all information and warnings in the documentation dealing with the products used together with the sensor (housings, transmitters, etc.).
- observing all safety regulations governing the equipment in which the sensor is installed.
- observing the prescribed inspection and maintenance intervals.
- correct care and maintenance of the unit, according to the instruction manual.
- observing the prescribed environmental and operational conditions, and permitted installation positions.

2.3 Safety instructions

- The InPro 5000 (i) sensors should be installed, operated, maintained only by personnel familiar with the sensor and who are qualified for such work.
- A defective sensor must neither be installed nor put into service.

- Only the maintenance work described in these operating instructions may be performed on the sensors.
- When changing faulty components, use only original spare parts obtainable from your METTLER TOLEDO supplier (see section 9).
- No modifications to the sensors and the accessories are allowed. The manufacturer/supplier accepts no responsibility for damage caused by unauthorised modifications. The risk is borne entirely by the user.

2.4 Environmental protection

Waste electrical products should not be disposed of with household waste. Please recycle where facilities exist. Check with your Local Authority or retailer for recycling advice.

3 Product description

3.1 General information

The **sensor InPro 5000 (i)** with integrated temperature probe is used for dissolved carbon dioxide measurement.

The sensor has been designed for accurate and precise measurement at low to medium carbon dioxide partial pressure.

The sensor is **sterilizable** and **compatible with CIP** (cleaning in place).

The sensor consists of an exchangeable pH electrode which is built in a steel shaft and fixed with a small **hexagonal screw**.

3.2 Principles of CO₂ measurement

Here is a short summary of the principle of the potentiometric measurement of CO₂ which this sensor is based (Severinghaus). The InPro 5000 (i) sensor employs a gas permeable silicone membrane which is tightly stretched around a special engineered flat pH membrane. The CO₂ gas from the sample or process diffuses across the membrane until its partial pressure equilibrates within the electrolyte film in front of the pH sensitive glass and the bulk of the internal electrolyte.

The dissolved CO₂ gas reacts with water to form bicarbonate and H⁺ ions:

The formed H⁺ ions lead to a pH change of the inner electrolyte which is measured with the integrated pH electrode. Thus, the CO₂ partial pressure can be calculated using the pH and the temperature, which is detected as well.

3.3 Scope of delivery

Each sensor is supplied together with:

- 1 bottle CO₂-electrolyte 25 ml
- 1 quality control certificate
- 1 inspection certificate 3.1
(in accordance with EN 10204.3/3.1)

Membrane body and cap sleeve are not mounted.

The interior body is protected against drying out with a wetting cap, filled with CO₂-electrolyte.


Please check completeness.

3.4 Equipment features

InPro 5000

InPro 5000 i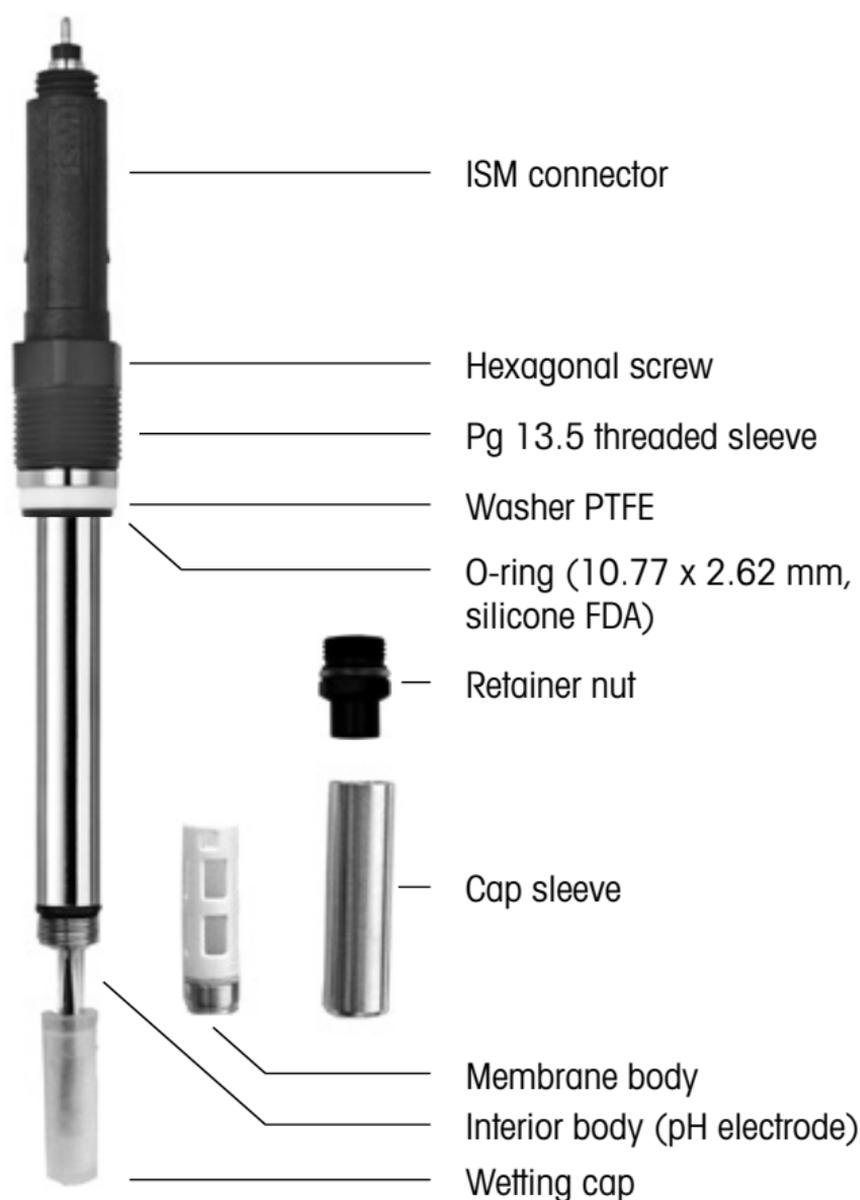


4 Preparation

4.1 Start-up

The interior body (pH electrode) is ready for calibration.

- On unpacking, check the pH electrode for mechanical damage. Report any damage immediately to your METTLER TOLEDO supplier.
- Remove the watering cap, cap sleeve, membrane body and briefly rinse the sensor with de-ionized water.
- After rinsing, the sensor should only be dapped dry with a tissue. Do not rub the pH sensitive glass, since this can lead to electrostatic charging and sluggish response times.
- Check inside the pH sensitive glass for the presence of any air bubbles and remove same by gently swinging in a vertical plane (shaking down as with a thermometer).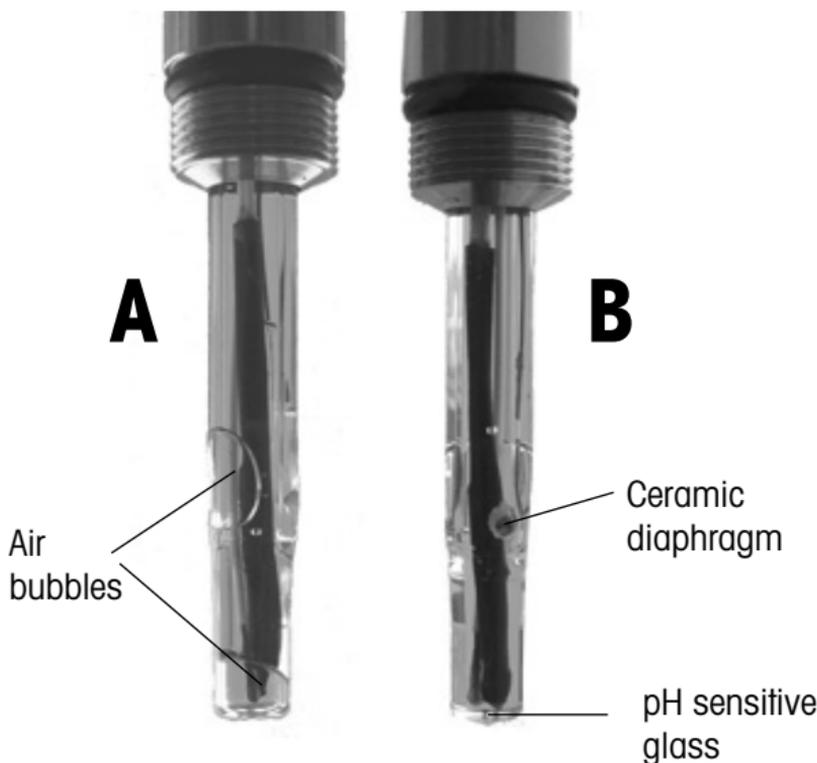

A: pH electrode with air bubbles will not work.


B: Correct pH electrode with removed bubbles.

Before using the sensor for the first time, make sure that the interior body (pH electrode) is well fixed into the steel body. The VP connector must not be rotated in the shaft. If so, the inner part has to be fixed by tightening the small **hexagonal screw** below the connector.

4.2 Calibration

Purpose of calibration: Each CO₂ sensor and Interior body (pH electrode) has its individual slope and zero point.


4.2.1 Connecting the InPro 5000 (i) to cable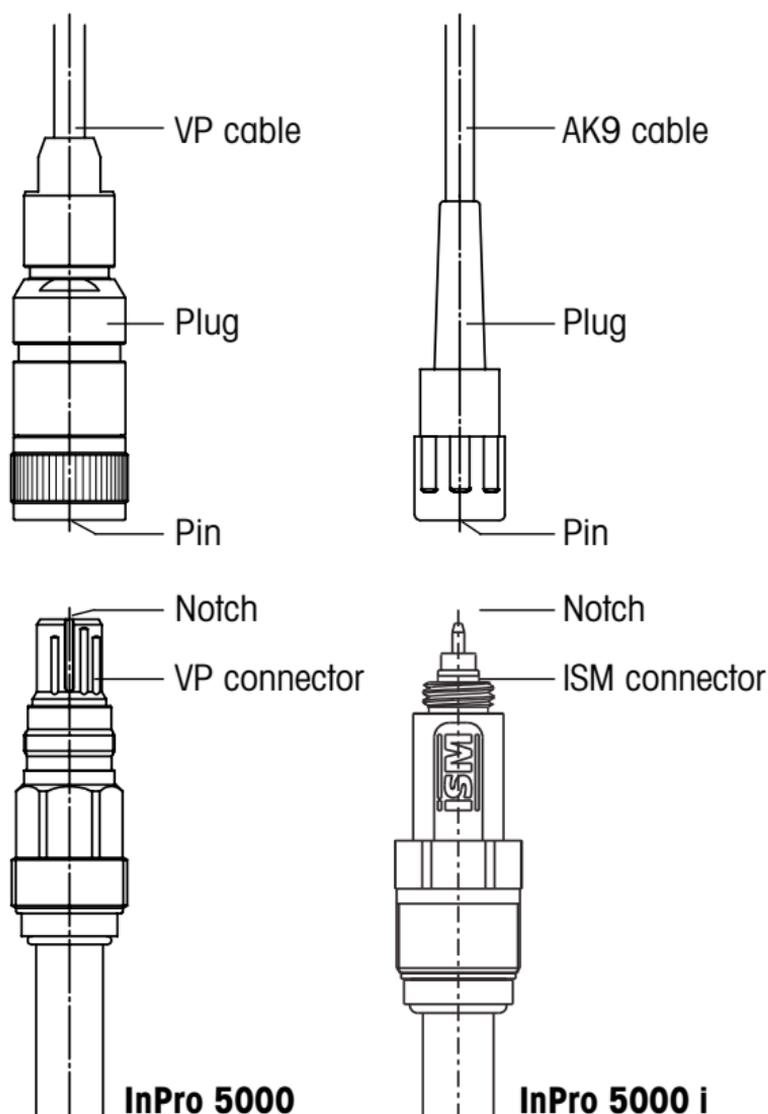

The sensor is connected to the transmitter via a VP or AK9 cable ensuring a secure connection under harsh industrial conditions. The robust watertight IP 68 connector housing guarantees maximum process safety.

To connect the VP cable to the sensor align the notch of the VP connector with the key in the plug. Then tightly screw the plug to fasten the two parts.


4.2.2 Connecting the cable to the transmitter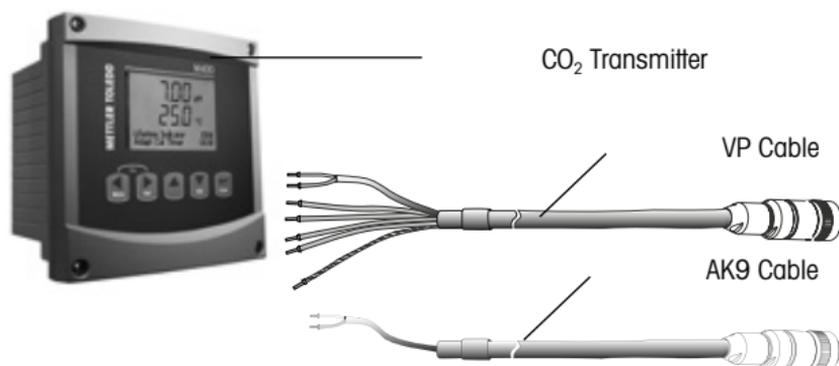

For connecting the sensor please refer to the manual of the transmitter.

4.2.3 Calibration of pH electrode

(Not needed for ISM digital InPro 5000 i)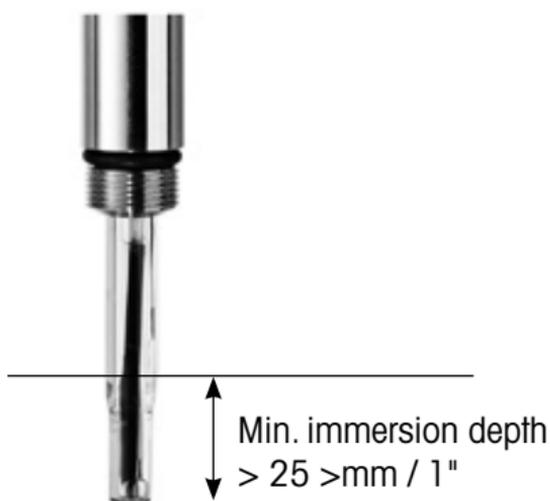


Prepare electrode as described in section 4.1 "Start-up". For the analog CO₂ sensor InPro 5000, we recommend a 2-point calibration in pH buffer pH 7.00 and pH 9.21. Before calibration, first remove wetting cap. Normally you start with buffer pH 7.00 for determining zero point and then buffer pH 9.21 for determining the slope. Please refer to instruction manual of the transmitter for further details.

Important settings : On the transmitter, select Pt1000 as temperature sensor for the analog InPro 5000. ISM digital InPro 5000 i does not require a temperature setting.

Important: Please do not leave the sensor in the calibration buffer for longer than 1 hour. Too long exposure may result in a high sensor drift in the process.

4.3 Mounting the membrane body

- After calibration of the pH electrode rinse the pH electrode with CO₂ electrolyte solution.
- Fill the new membrane body half way with CO₂ electrolyte.

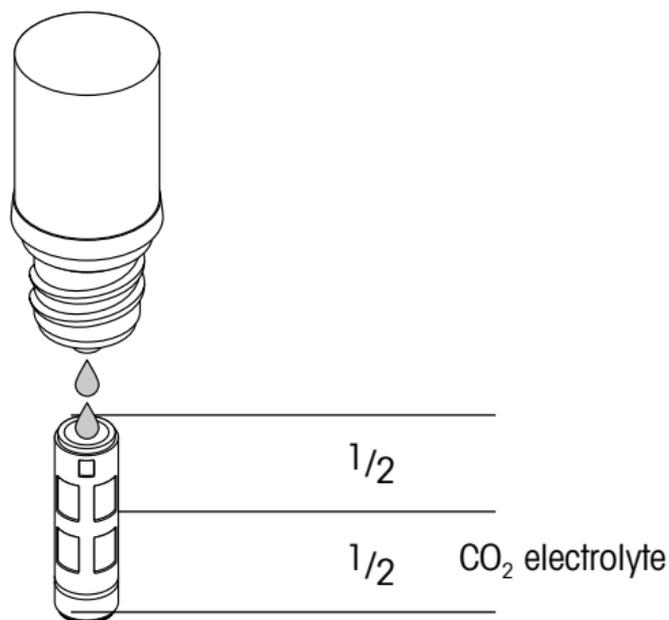Make sure that all air bubbles are removed from the membrane body. Air bubbles can be removed by carefully tapping on the membrane body.

- Slip the membrane body over the pH electrode while holding the sensor in a vertical position. The excess electrolyte will be displaced should be wiped dry with a paper tissue.

Important: No electrolyte, sample media or contamination may be present between the membrane body and the cap sleeve. Please check carefully.

- Carefully slip the cap sleeve over the fitted membrane body, holding the sensor in a vertical position and screw it tight. The cap sleeve must be clean and tight.

4.4 Mounting the sensor

Please read carefully the instruction manual of the particular housing (InFit761-... or InTrac797) for correct sensor installation.

The InPro 5000 (i) sensor can be mounted directly on a top plate of a small bioreactor or in a pipe with an inside Pg 13.5 thread and securely tightened via the Pg 13.5 threaded sleeve.

Remark: Mounting upside-down or horizontal is not possible due to internal pH electrode. **An angle of 15° or bigger to horizontal is needed.**

4.5 Sterilization

The installed sensor will be sterilized together with the bioreactor. The InPro5000 is in-situ sterilizable or in an autoclave. After cooling down but before inoculation a process calibration step, as described in chapter 5, is needed to achieve accurate measurements.

Remark: For best accuracy please **wait one hour after cooling down** before process calibration step according chapter 5.1.

5 Operation

5.1 Calibration of the CO₂ sensor

The InPro 5000 (i) sensor can be calibrated with CO₂ gas or a CO₂ / N₂ gas mixture. Example: If a CO₂ partial pressure of 150 mbar is expected, you can use a gas with 15 % CO₂/ 85 % N₂ in nitrogen to achieve best results. The calibration can be made in a gas or liquid phase. For calibration in a solution you have to wait until an equilibrium is reached. Depending on the pH value of the solution a large amount of CO₂ will react to build bicarbonate ions before building-up a constant partial pressure of CO₂.

Recommendations:

- **For small “benchtop” bioreactors** the sensor **can be calibrated by flushing the empty or liquid filled reactor with a CO₂ containing gas mixture.**
- **For larger bioreactors** the gas consumption would be too high and therefore **we recommend the calibration with the aid of the retractable housing InTrac 797.** Please operate the retractable housing according the instruction manual InTrac 797 to ensure sterile operation. Additionally the in- and outlet of the lower flushing chamber must be equipped with sterile filters to avoid any contamination with the calibration gas.

Calibration procedure:

- Sterilize CO₂ sensor InPro 5000 (i) in measuring position (membrane inside bioreactor) together with the bioreactor.
- Sterilize lower and upper flushing chambers of housing InTrac 797.
- After cooling down retract sensor in maintenance position (membrane in lower chamber).
- Flush lower chamber with sterile calibration gas. The pressure inside the flushing chamber must be known for proper calibration.
- After calibration close in- and outlet of the lower chamber and reinsert sensor in measuring position. Sensor is now ready for measurement.

6 Maintenance

6.1 Inspection

6.1.1 Visual inspection

To check the sensor, we recommend the following procedure:

- The contacts of the connector must be dry. Moisture, corrosion and dirt can lead to false readings.
- Check the cable for kinks, brittle areas or tears.
- Examine the membrane foil optically for signs of damage. The foil must be intact and clean. Dirty membranes should be wiped clean using a soft, moist tissue.

Note: An undulated membrane has no influence on the sensor performance, assuming the membrane is intact.

- The membrane body must be replaced if the sensor has a long response time, the reading is unstable or subject to excessive drift, and if the sensor cannot be calibrated with CO₂ gas or the membrane shows sign of mechanical damage.
- Check the pH sensitive glass for cracks. The diaphragm should show a white color. If necessary rinse with deionized water and clean carefully with soft paper tissue.

Attention! Do not use any cleaning agents containing alcohol. This could damage the sensor or lead to fault results.

Attention! The pH electrode is made of glass and therefore fragile.

6.1.2 Checking wiring and transmitter with the pH simulator

Necessary equipment for analog CO₂ sensor InPro 5000:

pH simulator 112 (p/n 59 906 431) and VP module (p/n 52 120 939).

This equipment enables the testing of the VP cable and correct wiring of the cable to the transmitter. Switch the VP simulator to Pt 1000 position. Additionally it is possible to check the linearity, temperature compensation and quality of the input circuit (input resistance and input current) of the CO₂ transmitter.

6.2 Changing the CO₂ electrolyte and membrane body

We recommend to use a new membrane body after each batch (see section 4.3). After each calibration of the pH electrode (section 4.2) new CO₂ electrolyte should be filled in the membrane body.

6.3 Maintenance of the interior body (pH electrode)

If the pH electrode shows a sluggish response time and / or insufficient sensitivity the pH sensitive glass can be reactivated using a HF solution (p/n 51 319 053). The pH sensitive glass part only is immersed into this solution for 5 minutes, is then immediately and thoroughly rinsed with deionized water and then stored in CO₂ Electrolyte up to the diaphragm for 12 hours. This procedure will make the electrode faster but will limit the life time hardly.

6.4 Replacement of the interior body (pH electrode)

Typically after 10 – 20 sterilization cycles the pH electrode should be replaced. Please refer to the instruction manual of the replacement interior body InPro 5000 (i).

6.5 ISM

DLI: Dynamic Lifetime Indicator

The DLI provides information about the remaining lifetime of the interior body. As long as the DLI is above zero days the system is within the specified accuracy after a calibration. If the DLI is zero the inner body needs to be replaced.

Contributing factors for ageing:

- temperature during measurement
- number of CIP cycles
- number of SIP cycles
- number of autoclavings

The DLI is calculated in the following way.

Continuously: With the above parameters an actual sensor stress is calculated. During the operation the sensor load is increased. The accumulated sensor load divided by the elapsed time is the basis of the calculation of the remaining lifetime.

TTM: Time to Maintenance

The TTM gives an indication when maintenance needs to be performed. Maintenance includes

1. Electrolyte exchange
2. Cleanig of the inner body
3. Replacement of the membrane body

ACT: Adaptive Calibration Timer

The ACT provides information as to when the next calibration is required to ensure measurements will remain within the specified accuracy. This calculation is based on the DLI information.

Calibration history

The last four calibrations and the factory calibration data are stored in the sensor memory. These data can be read out with a transmitter or with the iSense Asset Suite software.

The calibration history gives valuable information regarding the quality of the calibration and the ageing of the interior body.

7 Storage

For storage we recommend to fill the membrane body with fresh CO₂ electrolyte and rebuilt the sensor. The wetting cap, filled with CO₂-electrolyte, will prevent a drying-out of the membrane.

8 Product specification

8.1 Certificates

Each sensor is delivered with a set of 3.1 certificates (complying with EN 10204/3.1).

All wetted metal parts (sensor shaft, cap sleeve and membrane body) are identified with an engraved symbol corresponding to **the heat number** on the paper certificate delivered with the sensor.

Each wetted metal part (sensor shaft, cap sleeve and membrane body) is polished in order to get a surface roughness lower than 0.4 µm (16 µin). This represents a roughness grade number of N5 (according to ISO 1320:1992).

8.2 Specifications

InPro 5000 (i)	
Measurement principle	Potentiometric / Severinghaus
Working conditions	
Mechanical pressure resistance	max. 3 bar at 25 °C [max. 42 psi at 77 °F]
Pressure range measurement	0.2 – 2 bar absolute [3 – 30 psi absolute]
Temperature range (measurement)	0...60 °C [32...140 °F]
Sterilization	< 130 °C [266 °F] in situ or autoclave
Construction	
Temperature signal	InPro 5000 i: Digital InPro 5000: Pt 1000
Cable	InPro 5000 i: AK9 InPro 5000: VP
O-ring material	Viton®, Silicone (FDA approved)
Membrane material	PTFE-silicone (reinforced)
Wetted metal parts	Stainless steel DIN 1.4435 similar to AISI 316 L
Dimensions	
Sensor diameter	12 mm (0.47")
Immersion length	120, 220, 320 mm (4.7, 8.66, 12.6")
Performance	
Measuring range	10 – 1000 hPa CO ₂
Accuracy	± 10 %, 600 – 900 hPa ≥ ± 10 %, > 900 hPa
Detection limit	10 hPa CO ₂
Response time	90 % of final value < 120 sec (at 25 °C [77 °F] from air to CO ₂)

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All rights reserved