# METTLER TOLEDO DG115-SC

**Glass electrode**

**Glaselektrode**

**Electrode de verre**

**Electrodo de vidrio**

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## **Combined pH electrode**

with ground-glass sleeve junction for direct pH measurements and acid/base titrations in aqueous solutions – for media with low ion concentration.

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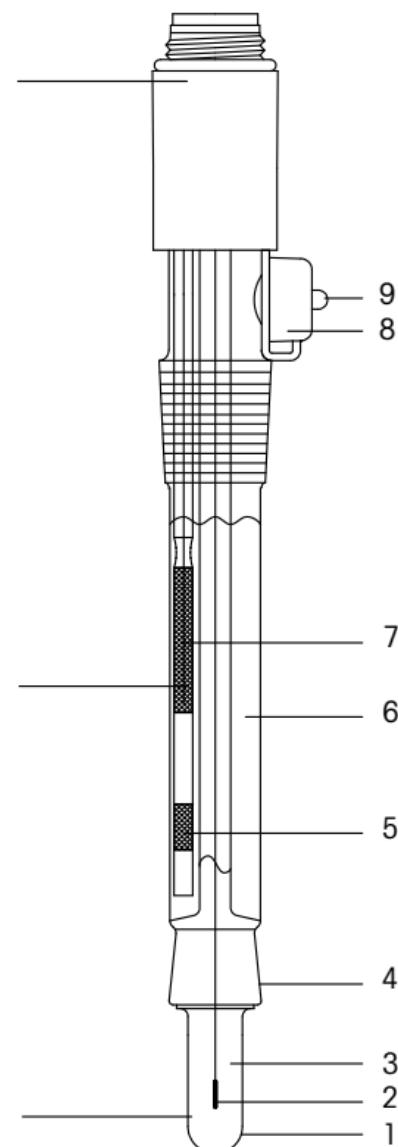
## **1. Start-up maintenance**

- 1.1 The internal buffer solution (3) must cover the whole inside of the glass membrane (1). Remove air bubbles by gently shaking in vertical direction.
- 1.2 To equalize pressure remove rubber cap (8) or perforate its nipple (9).
- 1.3 Refill the reference electrolyte (6) – (e.g. METTLER TOLEDO No. 51340045)– to approx. 1 cm below the aperture (8).
- 1.4 To activate the glass membrane place the electrode in a tenfold diluted reference electrolyte solution. Use a 0.5 mol/L KCl solution if no reference electrolyte is at hand.

## **2. Adjustment and calibration**

The parameters zero-point and slope are measured by adjustment. A new electrode should have parameter values according to 2.1–2.2. Instructions for the adjustment method are given in the titrator manual.

- 2.1 The zero point of the electrode (pH value at  $E = 0 \text{ mV}$ ) is approx. pH 7. The electrode potential, measured in a buffer solution of pH 7, must be within  $-30 \text{ mV}$  and  $+30 \text{ mV}$ .
- 2.2 The electrode slope (potential change per pH unit) at the pH range 4–9 must be greater than  $-55 \text{ mV/pH}$  (at  $25^\circ\text{C}$ ).
- 2.3 You should test the response time of the electrode. After 30 s the indicated electrode potential should not change more than 2 mV within the following 30 s.
- 2.4 If calibrating a used electrode the zero point should be within pH 6 and 8, the slope not less than  $-52 \text{ mV/pH}$ , and the potential change after 60 s should not exceed 3 mV/30 s.

## **3. Storage and lifetime**

- 3.1 Store the electrode in the reference electrolyte – the junction (4) has to be immersed – and close the aperture (perforated rubber cap is sufficient).
- 3.2 Never let the glass membrane dry out, otherwise it has to be reactivated and the KCl crystals in/on the junction have to be dissolved.
- 3.3 The lifetime of the electrode is 6 months to 3 years, depending on its use and treatment.

## **4. Maintenance instructions**

- 4.1 The measuring range extends from pH 0 to 14, the tolerable temperature range from 0 to 100 °C (32–212 °F).

- 4.2 The sample solution should never enter through the sleeve junction into the reference electrode. Therefore the level of the reference electrolyte must always be higher than the sample solution level.
- 4.3 No reaction should take place between the sample solution and the reference electrolyte: sulfides, bromides, iodides and cyanides react with  $\text{Ag}^+$  and some cations ( $\text{Ag}^+$ ,  $\text{Hg}^{2+}$ ,  $\text{Au}^{3+}$ ,  $\text{Pb}^{2+}$ ) react with  $\text{Cl}^-$  of the reference electrolyte and cause a blockage of the junction.
- 4.4 Solutions containing protein quickly block the junction, because protein precipitates on it.
- 4.5 Just wipe the electrode after rinsing, never rub it dry.
- 4.6 Do not let the cable connector plug corrode.
- 4.7 A dried out glass membrane can be reactivated by placing it for several hours in 0.1 mol/L HCl.

## 5. Cleaning procedures

- 5.1 Contamination of the reference electrolyte: Drain the reference electrolyte through aperture (8) and proceed according to 1.3 and 1.4.
- 5.2 Blocked junction by  $\text{AgCl}$ : place it overnight in concentrated ammonia, rinse and condition it for 1 hour in a buffer solution pH 4.
- 5.3 Blocked junction by sulfide precipitate: place it in a solution of thiourea (8% thiourea/1 mol/L HCl, e.g. METTLER TOLEDO No. 51340070) until the dark color has vanished.
- 5.4 Blocked junction by protein: place it at least for 1 hour in pepsin solution (5% pepsin/0.1 mol/L HCl, e.g. METTLER TOLEDO No. 51340068).
- 5.5 Other contaminations: place it for a few minutes in an ultrasonic bath with deion. water, ethanol or a mixed acid solution (e.g. aqua regia).

## 6. Regeneration

If the electrode does not meet the qualifications according to 2.4 anymore, try a corrosion of the glass membrane: place it 1–2 minutes in diluted HF solution (2% HF/5% HCl, e.g. METTLER TOLEDO No. 51340073), rinse thoroughly with deion. water and treat it as in 5.1.

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