**METTLER TOLEDO**

**DG114-SC**

**Glass electrode with Sensor Chip**

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**Plug & Play combined pH electrode**

with movable sleeve junction for direct pH measurements and titrations in aqueous solutions – for critical samples.

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# Care instructions

## Attention

Due to transportation and storage the movable sleeve ring of the sleeve junction may stick to the glass and thus stop the flow of electrolyte. This may cause a very instable or drifting signal and may increase the response time.

## Procedure

1. Close the filling aperture with the rubber cap.
2. Warm the movable sleeve ring with warm water and loosen it by hand by turning
3. Turn the electrode upside down and allow it to cool (loose sleeve junction).
4. Close the movable sleeve by a slight turning motion.
5. Check the fill level of the electrolyte and refill if necessary.
6. The electrode is ready to use.

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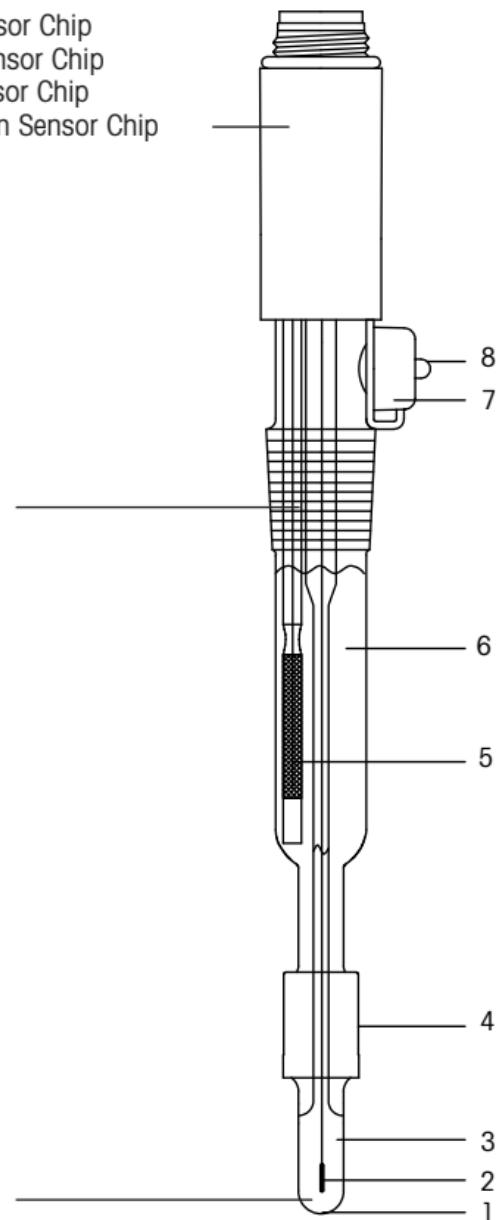
## **1. Start-up maintenance**

- 1.1 The internal buffer solution (3) must fill the entire inside of the glass membrane (1). Remove air bubbles by shaking the electrode gently.
- 1.2 To equalize pressure remove the rubber cap (7) or perforate its nipple (8).
- 1.3 Refill the reference electrolyte (6) – (e.g. ME-51340045) – just up to the aperture (7). Ensure that the silver chloride reference element (5) is always immersed in the reference electrolyte (6).
- 1.4 To activate the glass membrane place the electrode in a tenfold diluted reference electrolyte solution. Use a 0.5 mol/L KCl solution if no reference electrolyte is at hand.

## **2. Plug & Play performance**

The Sensor Chip which stores the sensor parameter is located in the screw cap of the sensor.

The sensor is recognized after connection to the sensor input of the titrator as well as after each restart of the system by reading the sensor parameter on the chip. A popup appears on the titrator terminal when the sensor was successfully recognized. This popup allows you to directly enter the sensor setup. You find further information concerning the functionality of the Plug & Play sensor.

## **3. Adjustment and calibration**

The parameters zero-point and slope are measured by calibration. A new electrode should have parameter values according to 3.1 – 3.2. Instructions for the calibration method are given in the titrator manual.

- 3.1 The zero-point of the electrode (pH-value at  $E = 0 \text{ mV}$ ) is approx. pH 7. The electrode potential, measured in a buffer solution of pH 7, must be between  $-30 \text{ mV}$  and  $+30 \text{ mV}$ .
- 3.2 The electrode slope (potential change per pH-unit) at the pH range 4–9 must be greater than  $-55 \text{ mV/pH}$  (at  $25^\circ \text{C}$ ).
- 3.3 You should test the response time of the electrode. After 30 s the indicated electrode potential should not change more than 2 mV within the following 30 s.
- 3.4 If testing a used electrode the zero-point should be between pH 6 and 8, the slope not less than  $-52 \text{ mV/pH}$ , and the potential change after 60 s should not exceed 3 mV/30 s.

#### 4. Storage and lifetime

- 4.1 Store the electrode in the reference electrolyte – the sleeve junction (4) has to be immersed – and close the aperture (perforated rubber cap is sufficient).
- 4.2 Never let the glass membrane dry out, otherwise it has to be reactivated and the KCl crystals in/on the sleeve junction have to be dissolved.
- 4.3 The lifetime of the electrode is 6 months to 3 years, depending on its use and treatment.

#### 5. Maintenance instructions

- 5.1 The measuring range extends from pH 0 to 14, the tolerable temperature range from 0 to 60 °C (32–140 °F).
- 5.2 The sample solution should never enter through the sleeve junction into the reference electrode. Therefore the level of the reference electrolyte must always be higher than the sample solution level.
- 5.3 Just blot the electrode after rinsing, never rub it dry.
- 5.4 Do not let the cable connector plug corrode
- 5.5 A dried out glass membrane can be reactivated by placing it for several hours in 0.1 mol/L HCl.

#### 6. Cleaning procedures

- 6.1 Contamination of the reference electrolyte: Drain the reference electrolyte through the aperture (7) and proceed according to 1.3 and 1.4.
- 6.2 Sleeve junction blocked by AgCl: place it overnight in concentrated ammonia, rinse and condition it for 1 hour in a buffer solution pH 4.
- 6.3 Sleeve junction blocked by sulfide precipitate: place it in a solution of thiourea (8% thiourea/1mol/L HCl, e.g. ME-51340070) until the dark color has vanished.
- 6.4 Sleeve junction blocked by protein: place it for at least 1 hour in pepsin solution (5% pepsin/0.1 mol/L HCl, e.g. ME-51340068)
- 6.5 Other contaminations: place it for a few minutes in an ultrasonic bath with deion. water, ethanol or a mixed acid solution (e.g. aqua regia).

#### 7. Regeneration

If the electrode does not meet the qualifications according to 3.4 anymore, try a corrosion of the glass membrane: place it 1–2 minutes in diluted HF-solution (2% HF/5% HCl, e.g. ME-51340073), rinse thoroughly with deion. water and treat it as in 6.1.

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